



DrosoFiller Replacement Teflon Plate Installation Instructions*

*Depending on the age of your filler you may want to replace old springs (cat#59-177) and add o-rings (cat#59-175)

- 1. Turn the DrosoFiller upside down exposing the bottom multi-hole steel plate.**
- 2. Using a long +/Phillips type screw driver carefully remove the screws, washers, springs, and o-rings - holding the plates together. We recommend placing these parts in a small bowl...do not lose any of these parts!**
- 3. Next remove the steel and Teflon plates. By design the fit of the steel plate can be tight over the posts. Slowly and gradually the plate will release.**
- 4. We recommend cleaning the bottom of the filler and steel plate before installing the new Teflon plate.**

Continued on back

5. Next place the new Teflon plate over the posts. The correct fit is, with your hand slide the plate back and forth. Notice the alignment of Teflon plate holes and holes on the bottom of the filler. If necessary flip the Teflon plate over for a better fit and alignment of holes.
6. Next reinstall the steel plate. Flat edge side facing the Teflon plate.
7. Next reinstall in this order - the o-ring, spring, washer, and screw.
8. Done! Call/e-mail FlyTabs with any questions.

flytabs@yahoo.com

805-748-5665