

# Bovine Growth Serum, US Origin

HYCLONE SERA

HyClone™ Bovine Growth Serum (BGS) was developed for cell culturists as an affordable and viable alternative to fetal bovine serum (FBS). Growth promotion studies demonstrate that BGS is an excellent replacement for FBS. BGS consists of high-quality bovine calf serum supplemented with chemically defined components including vitamins, amino acids, trace metals, and other small molecules to stimulate cell growth and proliferation.

BGS is filtered through three sequential 100 nm (0.1 µm) pore size-rated filters. Before dispensing, each lot of serum is pooled using true pool technology to ensure uniformity and consistency between bottles (Fig 1).

Key features of Bovine Growth Serum, US Origin

- Cost-efficient alternative to FBS
- Sourced in the USA
- Complete traceability back to original source
- Low in antibodies and high in growth factors
- Virus panel testing according to 9 CFR 113.53

## Product storage and handling

Sera should be stored at -10°C or lower. Once thawed, sera should be stored at 2°C to 8°C for up to six weeks in order to maintain quality. If the serum needs to be stored longer than six weeks after opening, it is recommended to aliquot the serum into convenient volumes and refreeze. Handle bottles that have been stored in freezer carefully. Avoid large temperature shifts and protect the serum from exposure to light. Refer to safety data sheet for any safety recommendations. Storage requirements are listed on the product label.

## Thawing

Remove serum from storage at -10°C or lower and place in a refrigerator overnight at 2°C to 8°C. Transfer the serum to a 37°C water bath, agitate periodically to mix the solutes concentrated at the bottom of the container. Do not hold the serum at 37°C any longer than necessary after thawing. Thawing serum in a bath above 40°C without mixing can denature the concentrated proteins in the bottom of the container and precipitates might form in the bottle. Thawing serum at higher temperatures is not recommended.



Fig 1. Bovine Growth Serum, US Origin.

Alternatively, bottles can be placed directly from storage at -10°C or lower into a 37°C water bath. Bottles should be agitated to enhance mixing and thawing. Turbidity and flocculent material might be present after thawing or after prolonged storage.

Experience indicates that regardless of the method used to thaw serum, it is critical that it is mixed during the thawing process to prevent the formation of gradients and subsequent precipitation. Because of differences in thawing rates of different components, serum will form a gradient if it is not mixed as it thaws. If serum is allowed to remain in such a gradient state, precipitation is likely to occur.

## General culture recommendations

Supplementation of classical media such as Dulbecco's Modified Eagle's Medium (DMEM) is recommended at a range between 5% and 10% BGS to support culture of a wide variety of cell lines and applications.

## Quality control testing

BGS is assayed for insulin, gamma globulin, alkaline phosphatase, lactate dehydrogenase, glutamic pyruvic transaminase (SGPT), glutamic oxaloacetic transaminase (SGOT), pH, total protein, albumin, blood urea nitrogen, creatinine, total bilirubin, sodium, potassium, calcium, chloride, phosphorus, osmolality, iron, total iron binding capacity (TIBC), percent saturation, glucose, and IgG. Assays are subject to change without notice.

Test specifications are listed in Table 1. Cell lines successfully cultured with BGS are listed in Table 2.

**Table 1.** Test specifications

Endotoxin (Limulus ameocyte lysate gel clot assay)	≤ 10 EU/mL
Hemoglobin (spectrophotometric)	≤ 10 mg/dL
Sterility testing (current USP)	
Bacteria and fungi	No growth
Virus testing (9 CFR 113.53)	
Fluorescent antibody	
Bluetongue	Not detected
Bovine adenovirus	Not detected
Bovine parvovirus	Not detected
Bovine respiratory syncytial virus	Not detected
Bovine viral diarrhea virus	Not detected
Rabies	Not detected
Reovirus	Not detected
Cytopathogenic agents (e.g., IBR)	Not detected
Hemadsorbing agents (e.g., PI3)	Not detected
Mycoplasma	
Large volume, direct culture	Not detected
Hoechst DNA stain	Not detected
Certificate of suitability	Included

**Table 2.** Cell lines successfully cultured with BGS

Cell Line	Description
COS-1/COS-7	African green monkey kidney fibroblasts
TT	Human epithelial thyroid medulla carcinoma cells
DMS53	Human lung carcinoma cells
HUVCC	Human umbilical vein epithelial cells
HeLa	Human epithelial adenocarcinoma cells
BHK	Baby hamster kidney fibroblasts
HEK293	Human kidney cells
Jurkat	Human thymoma cells
Hep G2	Human epithelial liver carcinoma cells
MRC5	Human lung fibroblasts
721	Mouse liver epithelial cells
HCT116/RKO	Human colorectal carcinoma cells
MCF-7	Human mammary gland adenocarcinoma cells
McCoy	Mouse fibroblasts
NT2	Human neuronal cells
PC-3	Human prostate carcinoma cells
WiDr	Human colorectal adenocarcinoma cells
BxPC-3	Human pancreatic adenocarcinoma cells
U-87	Human glioblastoma cells
CaSki	Human cervical carcinoma cells
T24	Human bladder carcinoma cells

## Related products

### ***HyClone classical media***

HyClone classical media are manufactured using ISO 9001- and ISO 13485-certified processes. All raw material components have passed strict quality control testing to ensure the appropriate level of quality. The classical media are hydrated using purified process water and have undergone 0.1 µm sterile filtration.

### ***HyClone phosphate buffered saline (PBS)***

Our PBS products are manufactured according to cGMP guidelines using ISO 9001- and ISO 13485-certified processes. The products have full traceability and documented origin of all formula ingredients.

### ***HyClone trypsin protease***

Our trypsin protease is derived from porcine pancreas and is gamma irradiated prior to hydration and filling. The product is formulated without calcium and magnesium.

## Ordering information

<b>Product</b>	<b>Size</b>	<b>Product code</b>
HyClone Bovine Growth Serum, US Origin	100 mL	SH30541.02
	500 mL	SH30541.03
	1000 mL	SH30541.04

Find certificates of analysis, safety data sheets, standard formulations, product inserts, and protocols at [www.cytiva.com/hyclonecerts](http://www.cytiva.com/hyclonecerts).

## **cytiva.com/hyclone**

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. HyClone is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](http://cytiva.com/contact)

CY14569-01Oct20-DF

