VORTEX-GENIE™ Series

VORTEX-GENIE™ 2

VORTEX-GENIE™ 2T

Digital VORTEX-GENIE™ 2

ABI Digital VORTEX-GENIE™ 2

VORTEX-GENIE™ Pulse

Instruction Manual

EN

ES

FR

DE

IT

Change History

#	Date	Version	Descriptions
1	2026.02.24	A	New release

Table of Contents

Introduction	2
Applicable Products	2
Safety Information	2
Package Contents	3
Intended Use	3
Installation	4
Specifications	5
Control Panel	
<i>VORTEX-GENIE 2</i>	6
<i>VORTEX-GENIE 2T</i>	6
<i>Digital VORTEX-GENIE 2 / ABI Digital VORTEX-GENIE 2</i>	7
<i>VORTEX-GENIE Pulse</i>	7
Applications	8
Operation Instruction	
<i>VORTEX-GENIE 2</i>	8
<i>VORTEX-GENIE 2T</i>	9
<i>Digital VORTEX-GENIE 2 / ABI Digital VORTEX-GENIE 2</i>	11
<i>VORTEX-GENIE Pulse</i>	13
Maintenance	15
Technical Support Information	15
Technical Data	16
Compliance	16

Introduction

This manual contains installation, operation, and maintenance instructions for OHAUS VORTEX-GENIE™ vortex mixers. Please read the manual completely before using.

Applicable Products

This instruction manual is applicable to the following products:

Product	Model Name
VORTEX-GENIE 2	e-V2 (SI-xxxx*)
VORTEX-GENIE 2T	e-V2T (SI-xxxx*)
Digital VORTEX-GENIE 2	e-V2D (SI-xxxx*)
ABI Digital VORTEX-GENIE 2	e-V2A (SI-xxxx*)
VORTEX-GENIE Pulse	e-V2P (SI-xxxx*)

* SI-xxxx represents the original SI model name.

Safety Information

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.

CAUTION For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.

ATTENTION For important information about the product. May lead to equipment damage if not avoided.

NOTE For useful information about the product.

Warning Symbols

General hazard

Electric shock hazard

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift unit by its attachment.

Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly.

DO NOT immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

DO NOT position the equipment such that it is difficult to disconnect the power cord during use. Always place the unit on a clean and sturdy work surface.

WARNING! The protection provided by the unit may be impaired if used with accessories not provided or recommended by the manufacturer or used in a manner not specified by the manufacturer.

WARNING! DO NOT use the unit in a hazardous atmosphere or with hazardous materials for which the unit was not designed.

CAUTION: To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplugging from the wall outlet.

Earth Ground – Protective Conductor Terminal

Alternating Current

Package Contents

- Vortex mixer
- Cup Head (not included in **ABI Digital VORTEX-GENIE 2**)

We recommend you retain the original packaging for 90 days in case you need to return the product for any reason to your distributor or Ohaus.

Intended Use

This instrument is intended for use in laboratories, pharmacies, schools, businesses and light industry. It must only be used for processing materials as described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use. If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

Installation

Upon receiving the unit, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred during transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the unit on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand the typical heat produced by the unit.

Do not position the equipment such that it is difficult to disconnect the power cord during use.

Always place the unit on a clean and sturdy work surface.

The unit is supplied with a power cord that should be plugged into a properly grounded outlet. If the supplied power cord does not meet your needs, please use an approved power cord that suits local codes and electric supply. Replacement of the cord must be made by a qualified electrician.

Specifications

Product	VORTEX-GENIE 2	VORTEX-GENIE 2T	Digital VORTEX-GENIE 2	ABI Digital VORTEX-GENIE 2	VORTEX-GENIE Pulse
Control	Analog		Digital		
Display	--		Red LED		
Working Mode	Touch, Hands Free				Hands Free, Pulse
Accessories Included	3-inch Platform, Pop-off Cup			Applied Biosystem's 96/384-Well Microplate Adaptor	3-inch Platform, Pop-off Cup
Speed Range	600 rpm - 3200 rpm 600 rpm - 2700 rpm (for 50Hz models)		500 rpm - 3000 rpm 500 rpm - 2850 rpm (for 50Hz models)		
Orbit	4 mm		3.5 mm		3 mm
Timer	--	1-60 seconds; 1-60 minutes	1-99 seconds; 1-99 minutes		1-99 minutes
Pulse - ON/OFF Time	--				0.1 second – 99 minutes
Power	120V, 0.65A, 60Hz 230V, 0.5A, 50Hz 100V, 1.0A, 50Hz		120V, 0.65A, 60Hz 230V, 0.5A, 50Hz 100V, 1.0A, 50/60Hz		
Dimensions (L x W x H)	165 x 122 x 165 mm / 6.5 x 4.8 x 6.5 inch				
Net Weight	4 kg / 8.8 lb.				

Control Panel

VORTEX-GENIE 2

A	Speed Dial
B	Speed Control Knob
C	Toggle Switch

VORTEX-GENIE 2T

A	Speed Dial
B	Timer Dial
C	Speed Control Knob
D	Timer Control Knob
E	Toggle Switch

Digital VORTEX-GENIE 2 / ABI Digital VORTEX-GENIE 2

A	Time Display
B	Speed Display
C	Mode / Timer Indicator
D	Mode Button
E	Timer Up / Down Buttons
F	Start / Stop Button
G	Speed Up / Down Buttons

VORTEX-GENIE Pulse

A	Time Display
B	Speed Display
C	Pulse On / Off Indicator
D	Pulse Button
E	Timer Up / Down Buttons
F	Start / Stop Button
G	Speed Up / Down Buttons

Applications

The 3-inch Platform accepts single or multiple test tubes, beakers, and various flasks. The removable cover, made of thermoplastic rubber, is highly impervious to chemicals. It is, however, recommended that spills be blotted quickly. Do not attempt to use the 3-inch Platform without the thermoplastic cover.

The Pop-off Cup can accommodate tubes of virtually any size or shape from micro-tubes to those up to 40mm in diameter and 200mm in length. In most instances, tubes without caps may be mixed with proper speed control. Tubes must be capped when potential aerosols may be hazardous.

The Microplate Adaptor for Applied Biosystems' Microplate Protocols securely holds Applied Biosystems' 96-well plates and 384-well plates.

Visit OHAUS.com for the complete accessory attachment options.

To install the Pop-off Cup, the 3-inch Platform, or any other accessory attachments, first align the "flat" part of the mixer shaft with the "flat" part of the attachment. **YOU MUST PRESS DOWN HARD WITH YOUR HAND UNTIL YOU HEAR THE ATTACHMENT SNAP INTO PLACE.** This will ensure that the attachment does not come off the mixer during operation.

Operation Instructions

VORTEX-GENIE 2

- 1.0 - Plug the cord into a properly grounded, three-pronged outlet.
- 2.0 - The VORTEX-GENIE 2 mixer has the following modes of operation.
- 2.1 - TOUCH MODE:

Place the toggle switch in the TOUCH position and set the speed control to any position between 3 and 10. A setting of 3 is least vigorous mixing and 10 is the most vigorous mixing. When you touch a vessel to the 3-inch Platform or Popoff Cup, the mixer will start automatically. Releasing the pressure on the Platform or Cup will stop the mixer.

2.2 - HANDS FREE MODE:

Place the toggle switch in the ON position and set the speed control to any position between 3 and 10. Hands Free mode allows for unattended operation that will continue indefinitely until you stop the mixer by placing the toggle switch into the OFF position. In this mode, you may use any of the accessory attachments available for the VORTEX-GENIE 2 mixer. **Review instruction No. 3.0 before attempting to change attachments.**

Note: The speed control switch should be set between positions 3 and 10 to enable the mixing action to start. After the action has started, you may, if desired, reduce the speed below position 3 for the duration of the run. Before re-starting the mixer, re-set the speed control to a setting higher than position 3.

- 3.0 - To change accessory mixing attachments, the following sequence is to be followed explicitly.
 - 3.1 - Turn the power OFF by placing the toggle switch into the OFF position.
 - 3.2 - Grasp the installed attachment and pull straight up until it releases from the shaft. A technique of lifting the attachment with your fingertips while using your knuckles as a fulcrum will give you the leverage to remove the installed attachment component smoothly.
 - 3.3 - Hold the part you wish to install so the "flat" part of the shaft is aligned with the "flat" of the receptacle in the attachment. Alternatively, you may place the attachment on the shaft and turn it until you feel the "flats" aligning.
 - 3.4 - Press attachment firmly downward until it "snaps" into place. The VORTEX-GENIE 2 mixer is now ready to use with the newly installed accessory attachment.
- 4.0 - With prolonged use, your VORTEX-GENIE 2 mixer may become warm to the touch. The motor will radiate a certain amount of heat that will dissipate by the cast zinc housing.

VORTEX-GENIE 2T

- 1.0 - Plug the cord into a properly grounded, three-pronged outlet.
- 2.0 - The VORTEX-GENIE 2T mixer has the following modes of operation.
 - 2.1 - "TT" (TIMED TOUCH - 1 to 60 seconds)

For procedures that require **short duration vortex** mixing from 1 to 60 seconds. Or, for procedures that require **accurate and repetitive vortex mixing sequences** of 1 to 60 seconds each.

In the "TT" operating mode, place the toggle switch in the TOUCH position and set the 1 to 60 second timer (gray knob with white reference line) to the desired mixing time on the gray numbered scale. Next, set the speed control (black pointer shaped knob) to any position between 3 and 10 on the black numbered scale before initiating the "TT" operating mode. A setting of 3 is the least vigorous mixing and 10 is the most vigorous mixing. When you touch a tube or other vessel to the 3-inch Platform or Pop-off Cup, the mixer will start and run for 1 - 60 seconds, depending on the timer setting. The mixer will automatically turn off at the end of the 1 – 60 second cycle. Releasing the pressure on the

Platform or Cup and then re-applying the pressure will initiate another 1 - 60 second cycle. The mixing cycles will be precisely timed, and you may initiate as many mixing cycles as you wish.

2.2 - "CT" (*CONTINUOUS TOUCH*)

The "CT" mode is used when you desire to operate the VORTEX-GENIE 2T so that vortex mixing starts and runs continuously, without interruption, when you touch the vessel to either the 3-inch Platform, or to the Pop-off Cup. To set the "CT" operating mode, place the toggle switch in the TOUCH position and turn the gray. Timer knob completely counterclockwise until it "clicks" into the continuous (NO TIMER) position. Set the speed control knob. (see 2.1 on previous page).

2.3 - "FT" (*HANDS FREE TIMED - 1 to 60 minutes*)

The "FT" mode allows for unattended operation that is timed from 1 to 60 minutes. In this mode, you may use any of the accessory attachments that are available for the VORTEX-GENIE 2T mixer. **Review instruction No. 3.0 before attempting to change attachments.**

To initiate the "FT" operating mode, place the toggle switch in the center (unmarked) position. With the desired accessory attachment installed and with the sample tubes or vessels in place, set the 1-to-60-minute timer (gray knob and gray numbered scale) to the desired run time. With the toggle switch in the center position, the VORTEX- GENIE 2T is OFF until the toggle switch is depressed to the START/ STOP position, held for a second and then released. This will start the mixing action of the "GENIE™". The mixer will run until the set time has elapsed. If for any reason you wish to stop the mixing action before the set time has elapsed, simply depress the toggle switch to the START/STOP position for a second time and release. The mixer will stop. Depressing the toggle switch a third time will initiate a completely new timed cycle.

2.4 - "FC" (*HANDS FREE CONTINUOUS*)

The "FC" mode allows for unattended operation that will continue indefinitely until you stop the run by depressing the toggle switch. In this mode, you may also use any of the accessory attachments available for the VORTEX-GENIE 2T mixer. To initiate the "FC" mode, turn the gray timer knob completely counterclockwise until it "clicks" into the continuous (NO TIMER) position. Next, depress the toggle switch to the START/STOP position for one second and then release. The VORTEX-GENIE 2T mixer will start and continue to run, unattended, until you once again depress the toggle switch to the START/STOP position and hold it for a second before releasing. If you wish to restart the mixer once again, simply depress the toggle switch for a second and release.

Note: The speed control switch should be set between positions 3 and 10 to enable the mixing action to start. After the action has started, you may, if desired, reduce the speed below position 3 for the duration of the run. Before re-starting the mixer, re-set the speed control to a setting higher than position 3.

3.0 - To change accessory mixing attachments, the following sequence is to be followed explicitly.

- 3.1 - Turn the power OFF by first placing the toggle switch into the TOUCH position momentarily and then moving the toggle switch to the center (unmarked) position.
- 3.2 - Grasp the installed attachment and pull straight up until it releases from the shaft. A technique of lifting the attachment with your fingertips while using your knuckles as a fulcrum will give you the leverage to remove the installed component smoothly.
- 3.3 - Hold the part you wish to install so the "flat" part of the shaft is aligned with the "flat" of the receptacle in the attachment. Alternatively, you may place the attachment on the shaft and turn it until you feel the "flats" aligning.
- 3.4 - Press attachment firmly downward until it "snaps" into place. The VORTEX-GENIE 2T is now ready to use with the newly installed accessory attachment. Follow instructions 1.0 through 2.4 for proper use of the mixer.
- 4.0 - With prolonged use, your VORTEX-GENIE 2T mixer may become warm to the touch. The motor will radiate a certain amount of heat that will dissipate by the cast zinc housing.

Digital VORTEX-GENIE 2 / ABI Digital VORTEX-GENIE 2

1.0 - Plug the line cord into a properly grounded electrical outlet. Illuminated numbers will appear in the TIME and SPEED windows indicating that the unit is ON and ready for use.

2.0 - The DIGITAL VORTEX-GENIE 2 mixer has the following modes of operation:

2.1 - "TT" (*TIMED TOUCH - 1 to 99 seconds*)

For procedures that require **short duration vortex** mixing from 1 to 99 seconds. Or, for procedures that require accurate and repetitive vortex mixing sequences of 1 to 99 seconds each.

To initiate the "TT" mode, press the **MODE** button until **SEC** is illuminated beneath the word **TOUCH**. Press the UP/ DOWN arrow buttons beneath the **TIME** window to set the desired mixing time. Next, set the speed control by pressing the UP/DOWN arrow buttons beneath the **SPEED** window.

When you touch a tube or other vessel to the 3-inch Platform or Pop-off Cup, the mixer will start and run for 1 to 99 seconds, depending on the **TIME** setting. The mixer will automatically turn off at the end of the 1 to 99 second cycle. Releasing the pressure on the 3-inch Platform or Pop-off Cup and then re-applying the pressure will initiate another 1 to 99 second cycle. The mixing cycles will be precisely timed and you may initiate as many mixing cycles as you wish.

2.2 - "CT" (*CONTINUOUS TOUCH*)

The "CT" mode allows for operation that will continue indefinitely until you release pressure on the 3-inch Platform or Pop-off Cup.

To initiate the "CT" mode, press the **MODE** button until **SEC** is illuminated beneath the word **TOUCH**. Next, press and hold the DOWN arrow button beneath the **TIME** window until the illuminated display reaches "0". Upon releasing the button, the display will have two dash marks (--), indicating that no time limit is set. Set the speed by pressing the UP/DOWN arrow buttons beneath the **SPEED** window. Apply pressure to the 3-inch Platform or Pop-off Cup to initiate mixer. Release pressure to stop mixer.

2.3 - "FT" (*HANDS FREE TIMED - 1 to 99 minutes*)

The "FT" mode allows for unattended operation that is timed from 1 to 99 minutes. In this mode, you may use any of the accessory attachments that are available for the DIGITAL VORTEX-GENIE 2 mixer. **Review instruction 3.0 before attempting to change attachments.**

To initiate the "FT" mode, press the **MODE** button until **MIN** is illuminated above the START/STOP text. Next, press the UP/DOWN arrow buttons beneath the **TIME** window to set the desired mixing time. Next, set the speed control by pressing the UP/ DOWN arrow buttons beneath the **SPEED** window. Press the START/STOP button to begin the mixing action. The mixer will run until the set time has elapsed. To stop the cycle before the set time has elapsed, press the START/STOP button. To initiate a new timed cycle, press the START/STOP button again.

2.4 - "FC" (*HANDS FREE CONTINUOUS*)

The "FC" mode allows for unattended operation that will continue indefinitely until you stop the cycle by pressing the START/STOP button. In this mode, you may also use any of the accessory attachments available for the DIGITAL VORTEX-GENIE 2 mixer.

To initiate the "FC" mode, press the **MODE** button until **MIN** is illuminated above the START/STOP text. Next, press the DOWN arrow beneath the **TIME** window until the illuminated display reaches "0". Upon releasing the button, the display will have two dash marks (--), indicating that no time limit is set. Next, set the speed control by pressing the UP/DOWN arrow buttons beneath the **SPEED** window. Upon pressing the **START/STOP** button, the mixer will start and continue to run, unattended, until you once again press the **START/STOP** button. To restart the cycle, press the **START/ STOP** button again.

2.5 - "SLEEP"

The mixer will automatically go into a "SLEEP" mode if the mixer is plugged in but not used for 10 minutes. The **TIME** and **SPEED** illuminated displays will go blank and the **SEC** or **MIN** illuminated display will blink. Press the **START/STOP** button to "WAKE UP" the mixer.

Note: The speed control can be adjusted up or down during the mixing cycle, however, the TIME can only be adjusted before the cycle begins.

3.0 - To change accessory mixing attachments, the following sequence is to be followed explicitly.

- 3.1 - Grasp the installed attachment and pull straight up until it releases from the shaft. A technique of lifting the attachment with your fingertips while using your knuckles as a fulcrum will give you the leverage to remove the installed component smoothly.
- 3.2 - Hold the part you wish to install so the "flat" part of the shaft is aligned with the "flat" of the receptacle in the attachment. Alternatively, you may place the attachment on the shaft and turn it until you feel the "flats" aligning.
- 3.3 - Press attachment firmly downward until it "snaps" into place. The DIGITAL VORTEX-GENIE 2 is now ready to use with the newly installed accessory attachment. Follow instructions 1.0 through 2.5 for proper use of the mixer.
- 4.0 - With prolonged use, your DIGITAL VORTEX-GENIE 2 mixer may become warm to the touch. The motor will radiate a certain amount of heat that will dissipate by the cast zinc housing.

VORTEX-GENIE Pulse

- 1.0 - Plug the line cord into a properly grounded electrical outlet. Illuminated numbers will appear in the **TIME** and **SPEED** windows indicating that the unit is ON and ready for use.
- 2.0 - The VORTEX-GENIE PULSE mixer has the following features:
 - 2.1 - TIMED VORTEXING
The VORTEX-GENIE PULSE may be used as a normal (non-pulsing) vortex mixer. Press the PULSE button so that pulse is "OFF". The left display is set time (in minutes) till automatic stopping of the unit. The right display is set speed.

Note: When the unit is running, actual speed will be displayed which can vary based on load. For heavier loads on unit, desired setting speed may not be achievable.

"UP" represents the actual maximum speed (achievable under the given load conditions at that time). To activate the "UP" speed setting, press the UP arrow button beneath the **SPEED** window until maximum speed is displayed. Release the UP arrow button and press it again. "UP" will appear in the display window. Press the START/STOP button to initiate a mixing cycle.

2.2 - CONTINUOUS VORTEXING

Set TIME to "0" for continuous (non-timed) operation. Display will read "--"


The unit will run until the START/STOP button is pressed.

Set the speed by pressing the UP/DOWN arrow buttons beneath the SPEED window.

2.3 - PULSED VORTEXING

The PULSED VORTEXING will work in CONTINUOUS or TIMED modes. To initiate the PULSE “ON” mode, press the **PULSE** button until **ON** is illuminated beneath the word **PULSE**. The TIME and SPEED displays will begin flashing. When display is flashing, you are setting the Pulse “RUN” time (left display) and the Pulse “STOP” time (right display).

The TIME setting can range from 1/10th of a second to 99 minutes. When the rightmost decimal point is displayed, the setting is in seconds. When the rightmost decimal point is **not** displayed, the setting is in minutes. A setting of 0-1 second is adjusted in 1/10th of a second increments; 1-60 seconds are adjusted in 1 second increments; 1-10 minutes are adjusted in 5 minute (30 second) increments; 10-99 minutes are adjusted in 1 minute increments. (See Figure 1 for display readouts and corresponding time settings).

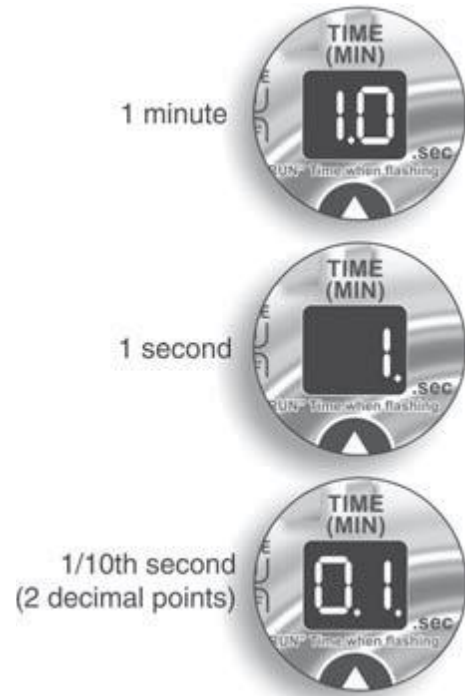Set the desired PULSE “RUN” time. Next set the desired PULSE “STOP” time. Wait for the TIME and SPEED displays to stop blinking.

It is recommended to set the **SPEED** to “UP” when using the PULSE mode with “RUN” times of less than 5 seconds (all speeds will work, however time/speed regulation is best at the “UP” setting).

Note: When using the PULSE mode with “RUN” times of less than 5 seconds, SET speed will be displayed when the unit is running, instead of actual speed (Minimum of 5 seconds is required to display a stable speed).

2.4 - “SLEEP”

The mixer will automatically go into a “SLEEP” mode if the mixer is plugged in but not used for 10 minutes. The **TIME** and **SPEED** illuminated displays will go blank and the PULSE **ON/OFF** illuminated display will blink. Press the **START/STOP** button to “WAKE UP” the mixer.

Note: The speed control can be adjusted up or down during the mixing cycle, however, the TIME can only be adjusted before the cycle begins.

3.0 - To change accessory mixing attachments, the following sequence is to be followed explicitly.

- 3.1 - Turn the power OFF by unplugging the unit from the outlet.
- 3.2 - Grasp the installed attachment and pull straight up until it releases from the shaft. A technique of lifting the attachment with your fingertips while using your knuckles as a fulcrum will give you the leverage to remove the installed component smoothly.
- 3.3 - Hold the attachment you wish to install so the "flat" part of the shaft is aligned with the "flat" of the receptacle in the attachment component. Alternatively, you may place the attachment component on the shaft and turn it until you feel the "flats" aligning.
- 3.4 - Press attachment component firmly downward until it "snaps" into place. The VORTEX-GENIE PULSE mixer is now ready to use with the newly installed accessory attachment. Plug the unit back into the electrical outlet. Follow instructions 1.0 through 2.4 for proper use of the mixer.
- 4.0 - With prolonged use, your VORTEX-GENIE PULSE mixer may become warm to the touch. The motor will radiate a certain amount of heat that will be dissipated by the cast zinc housing.

Maintenance

The OHAUS VORTEX-GENIE mixer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. DO NOT use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your Ohaus representative.

Technical Support Information

For service assistance or technical support, please visit our website at www.ohaus.com to locate and contact the Ohaus office nearest you.

Serial Number: _____

Date of Purchase: _____

Supplier: _____

Technical Data

Equipment Rating

- Indoor use only
- Altitude: 0 to 2000 m
- Environmental: 0°C to 38°C (32°F to 100°F), max 95% relative humidity, non-condensing
- Overvoltage category (Installation category): II

Compliance

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the applicable harmonized standards of EU Directives 2011/65/EU (RoHS), (EU)2015/863, 2014/30/EU (EMC) and 2014/35/EU (LVD). The EU Declaration of Conformity is available online at www.ohaus.com/ce .
	This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .
	EN 61326-1
	CAN/CSA-C22.2 No. 61010-1-12, CAN/CSA-C22.2 No. 61010-2-051 UL 61010-1, UL 61010-2-051

Global Notice

Warning: This is a Class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.

ISED Canada Compliance Statement:

CAN ICES-003(A) / NMB-003(A) FCC Notice

FCC Compliance Statement:

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

ISO 9001 Registration

The management system governing the production of this product is ISO 9001 certified.

FCC Supplier Declaration of Conformity

Unintentional Radiator per 47CFR Part B

Trade Name: OHAUS CORPORATION

Model: VORTEX-GENIE 2 e-V2..., VORTEX-GENIE 2T e-V2T..., Digital VORTEX-GENIE 2 e-V2D..., ABI Digital VORTEX-GENIE 2 e-V2A..., VORTEX-GENIE Pulse e-V2P...

Party issuing Supplier's Declaration of Conformity

Troemner LLC

201 Wolf Drive

Thorofare, NJ, 08086

USA

Phone: +1 856 686 1600

Responsible Party – U.S. Contact Information

Ohaus Corporation

8 Campus Drive, Ste. 105

Parsippany, NJ 07054

United States

Phone: +1 973 377 9000

Web: www.ohaus.com

LIMITED WARRANTY

OHAUS products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period OHAUS will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to OHAUS.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than OHAUS. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by OHAUS Corporation. OHAUS Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact OHAUS or your local OHAUS dealer for further details.

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P/N: 31018505 A © 2026 Ohaus Corporation, all rights reserved