

# F301.5 and FX301.5 Fixed-Angle Rotors

For Use in the Beckman Coulter Microfuge 22R (F301.5), Allegra X-22 Series (FX301.5), and Allegra X-30 Series (FX301.5) Centrifuges









#### F301.5 and FX301.5 Fixed-Angle Rotors

MMR-TB-002DB (September 2018)

© 2018 Beckman Coulter, Inc.

Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks mentioned herein are trademarks or registered trademarks of Beckman Coulter, Inc. in the United States and other countries.

All other trademarks, service marks, products, or services are trademarks or registered trademarks of their respective holders.

#### **Contact Us**

If you have any questions, contact our Customer Support Center.

- Worldwide, find us via our website at www.beckman.com/support/technical
- In the USA and Canada, call us at 1-800-369-0333.
- Outside of the USA and Canada, contact your local Beckman Coulter Representative.

Find us on the World Wide Web at: www.beckman.com

#### EC REP

Beckman Coulter Eurocenter S.A. 22, rue Juste-Olivier Case Postale 1044 CH - 1260 Nyon 1, Switzerland Tel: +41 (0) 22 365 36 11

Glossary of Symbols is available at beckman.com/techdocs (PN C24689).

**Original Instructions** 

## **Revision History**

This document applies to the latest and higher versions. When a subsequent version affects the information in this document, a new issue will be released to the Beckman Coulter website. For labeling updates, go to www.beckman.com/techdocs and download the latest version of the manual or system help for your instrument.

#### Issue DB 09/18

Changes were made to: Description; Care and Maintenance; Run Speeds.

**Note:** Changes that are part of the most recent revision are indicated in text by a bar in the margin of the amended page.

MMR-TB-002DB iii

IV MMR-TB-002DB

## Safety Notice

Read all product manuals and consult with Beckman Coulter-trained personnel before attempting to use this equipment. Do not attempt to perform any procedure before carefully reading all instructions. Always follow product labeling and manufacturer's recommendations. If in doubt as to how to proceed in any situation, contact your Beckman Coulter Representative.



This safety notice summarizes information basic to the safe use of the rotors described in this manual. The international symbol displayed to the left is a reminder to the user that all safety instructions should be read and understood before operation or maintenance of this equipment is attempted. When you see the symbol on other pages of this publication, pay special attention to the safety information presented. Observance of safety precautions will also help to avoid actions that could damage or adversely affect the performance of the rotor.

### Alerts for Warning, Caution, and Note



WARNING indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury.



CAUTION indicates a potentially hazardous situation, which, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices.

**NOTE** NOTE is used to call attention to notable information that should be followed during installation, use, or servicing of this equipment.

## Safety Information for the F301.5 and FX301.5 Rotors

Handle body fluids with care because they can transmit disease. No known test offers complete assurance that such fluids are free of micro-organisms. Some of the most virulent — Hepatitis (B and C) viruses, HIV (I–V), atypical mycobacteria, and certain systemic fungi — further emphasize the need for aerosol protection. Handle other infectious samples according to good laboratory procedures and methods to prevent spread of disease. Because spills may generate aerosols, observe proper safety precautions for aerosol containment. Do not run toxic, pathogenic, or radioactive materials in this rotor without taking appropriate safety precautions. Biosafe containment should be used when Risk Group II materials (as identified in the World Health Organization Laboratory Biosafety Manual) are handled; materials of a higher group require more than one level of protection.

MMR-TB-002DB V

The rotor and accessories are not designed for use with materials capable of developing flammable or explosive vapors. Do not centrifuge such materials in nor handle or store them near the centrifuge.

These rotors were developed, manufactured, and tested for safety and reliability as part of a Beckman Coulter centrifuge/rotor system. Their safety or reliability cannot be assured if used in a non-Beckman Coulter centrifuge, in a Beckman Coulter centrifuge not specified for use with the rotors, or in a Beckman Coulter centrifuge that has been modified without Beckman Coulter's approval.

Although rotor components and accessories made by other manufacturers may fit in the F301.5 and FX301.5 rotors, their safety in these rotors cannot be ascertained by Beckman Coulter. Use of other manufacturers' components or accessories in the F301.5 or FX301.5 rotors may void the rotor warranty and should be prohibited by your laboratory safety officer. Only the components and accessories listed in this publication should be used in these rotors.

Make sure that filled containers are loaded symmetrically into the rotor and that opposing tubes are filled to the same level with liquid of the same density.

If disassembly reveals evidence of leakage, you should assume that some fluid escaped the rotor. Apply appropriate decontamination procedures to the centrifuge and accessories.

Never exceed the maximum rated speed of the rotor and labware in use. Refer to the section on *Run Speeds*.

Do not use sharp tools on the rotor that could cause scratches in the rotor surface. Corrosion begins in scratches and may open fissures in the rotor with continued use.

Vİ MMR-TB-002DB

## Contents

```
Revision History, iii
Safety Notice, v
Alerts for Warning, Caution, and Note, v
Safety Information for the F301.5 and FX301.5 Rotors, {\bf v}
F301.5 and FX301.5 Fixed-Angle Rotors, 1
Specifications for the F301.5 Rotor, 1
Specifications for the FX301.5 Rotor, 2
Description, 3
Preparation and Use, 3
         Prerun Safety Checks, 3
Rotor Preparation, 4
Operation, 5
Removal and Sample Recovery, 5
         Tubes, 6
Run Speeds, 7
Care and Maintenance, 8
         Maintenance, 8
         Cleaning, 9
         Decontamination, 10
         Sterilization and Disinfection, 10
         Storage, 10
Returning a Rotor, 11
Supply List, 11
         Replacement Rotor Parts, 12
         Other, 12
Beckman Coulter, Inc.
Benchtop Rotor Warranty
```


## Tables

1 Available Tubes for the F301.5 and FX301.5 Rotors, 6

## F301.5 and FX301.5 Fixed-Angle Rotors

## **Specifications for the F301.5 Rotor**

**NOTE** The F301.5 rotor is used in the Microfuge 22R centrifuge only.

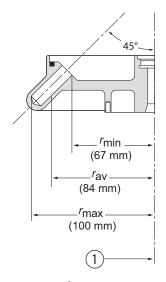

1. Axis of Rotation

Maximum speed
Density rating at maximum speed 1.2 g/mL
Critical speed range <sup>a</sup>
Relative Centrifugal Field <sup>b</sup> at maximum speed at $r_{\rm max}$ (100 min)
k factor at maximum speed
Conditions requiring speed reductions see Run Speeds
Maximum imbalance of opposing loads 6 grams
Number of tube cavities
Available bottles and tubes see Table 1
Nominal tube dimensions
Nominal tube capacity (largest tube)
Nominal rotor capacity
Approximate acceleration time to maximum speed (fully loaded)
Approximate deceleration time from maximum speed (fully loaded)
Weight of fully loaded rotor 1.67 kg (3.68 lb)
Rotor and lid material aluminum


a. The critical speed range is the range of speeds over which the rotor shifts so as to rotate about its center of mass. Passing through the critical speed range is characterized by some vibration.

MMR-TB-002DB 1-1

b. Relative Centrifugal Field (RCF) is the ratio of the centrifugal acceleration at a specified radius and speed  $(r\omega^2)$  to the standard acceleration of gravity (g) according to the following formula: RCF =  $r\omega^2/g$  — where r is the radius in millimeters,  $\omega$  is the angular velocity in radians per second  $(2 \pi \text{ RPM}/60)$ , and g is the standard acceleration of gravity (9807 mm/s<sup>2</sup>). After substitution: RCF =  $1.12r (\text{RPM}/1000)^2$ 

## **Specifications for the FX301.5 Rotor**

**NOTE** The FX301.5 rotor is used in the Allegra X-22 and Allegra X-30 Series centrifuges only.

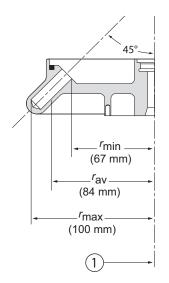

#### 1. Axis of Rotation


Maximum speed				
Allegra X-30R centrifuge				
Density rating at maximum speed 1.2 g/mL				
Critical speed range <sup>a</sup> 600 to 800 rpm				
Relative Centrifugal Field <sup>b</sup> at maximum speed at $r_{\rm max}$ (100 min) Allegra X-30R centrifuge 28,672 × $g$ Allegra X-22R centrifuge 21,952 × $g$ Allegra X-30 and Allegra X-22 centrifuges 19,515 × $g$				
k factor at maximum speed				
Conditions requiring speed reductions see Run Speeds				
Maximum imbalance of opposing loads 6 grams				
Number of tube cavities				
Available bottles and tubes see Table 1				
Nominal tube dimensions				
Nominal tube capacity (largest tube)				
Nominal rotor capacity				
Approximate acceleration time to maximum speed (fully loaded)				
Approximate deceleration time from maximum speed (fully loaded)				
Weight of fully loaded rotor				
Rotor and lid material aluminum				

a. The critical speed range is the range of speeds over which the rotor shifts so as to rotate about its center of mass. Passing through the critical speed range is characterized by some vibration.

1-2 MMR-TB-002DB

b. Relative Centrifugal Field (RCF) is the ratio of the centrifugal acceleration at a specified radius and speed  $(r\omega^2)$  to the standard acceleration of gravity (g) according to the following formula: RCF =  $r\omega^2/g$  — where r is the radius in millimeters,  $\omega$  is the angular velocity in radians per second (2  $\pi$  RPM /60), and g is the standard acceleration of gravity (9807 mm/s<sup>2</sup>). After substitution: RCF = 1.12r (RPM/1000)<sup>2</sup>

## **Description**

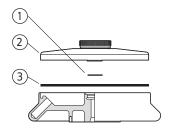

- 1. Lid O-ring (368896)
- 2. Lid (368895)
- **3.** O-ring (368991)

These Beckman Coulter rotors have been manufactured in an ISO 9001 or 13485 facility for use with the specified centrifuges.

The F301.5 and FX301.5 are fixed-angle rotors with a tube angle of 45 degrees from the axis of rotation. These rotors can centrifuge up to thirty 1.5 to 2.2-mL reaction-vials, as well as 250 to 400- $\mu$ L vials and 500 to 750- $\mu$ L tubes. These rotors develop centrifugal forces that can efficiently pellet subcellular organelles, viruses, bacteria, mitochondria, chloroplasts, or algae.

Although the rotors are similar in form and function, they are not interchangeable; the F301.5 rotor cannot be used in Allegra X-22 Series or Allegra X-30 series centrifuges, and the FX301.5 rotor cannot be used in the Microfuge 22R.

Refer to the Warranty at the back of this manual for warranty information

## **Preparation and Use**

Specific information about the F301.5 and FX301.5 rotors is given here. Information about the centrifuges is contained in the respective centrifuge manuals, which should be used together with this manual for complete centrifuge, rotor, and accessory operation.

**NOTE** Although rotor components and accessories made by other manufacturers may fit in the F301.5 or FX301.5 rotor, their safety in these rotors cannot be ascertained by Beckman Coulter. Use of other manufacturers' components or accessories in these rotors may void the rotor warranty and should be prohibited by your laboratory safety officer. Only the components and accessories listed in this publication should be used in these rotors.

## **Prerun Safety Checks**



Read the Safety Notice section at the front of this manual before using the rotor.

Make sure that the rotor, lid, and all tubes and accessories are clean and show no signs of corrosion or cracking.

- **2** If fluid containment is required, use capped tubes.
  - All containers carrying physiological fluids should be capped, and not overfilled, to prevent leakage.
  - Verify that the tubes being used are listed in Table 1.
- 3 Check the chemical compatibilities of all materials used (refer to *Chemical Resistances*, publication IN-175).

## **Rotor Preparation**

For runs at other than room temperature, refrigerate or warm the rotor before the run for fast equilibration.

- 1 Be sure that the metal threads in the rotor are clean and lightly but evenly coated with Spinkote lubricant (306812).
- **2** Load the filled containers symmetrically into the rotor (see *Tubes*, page 6 for tube information).
  - If fewer than 30 tubes are being run, they must be arranged symmetrically in the rotor.
  - Opposing tubes must be filled to the same level with liquid of the same density and weigh approximately the same (within 6 grams).
- **3** Put the lid in place and tighten it by hand.

1-4 MMR-TB-002DB

### **Operation**

Precool the rotor in the centrifuge or in a refrigerator before use—especially before short runs—to ensure that the rotor reaches the set temperature. A suggested precooling cycle is a minimum of 30 minutes at 2000 rpm with the temperature set to 0°C.

**NOTE** Temperatures may vary slightly between centrifuges. If sample temperature is crucial, test temperature settings on your instrument using water samples. For runs at other than room temperature, refrigerate or warm the rotor beforehand for fast equilibration.

Ensure that the rotor tie-down screw (365806 for the F301.5, 361367 for the FX301.5) is in good condition and the threads are free of foreign matter.



Do not drop the rotor onto the drive shaft. The shaft can be bent or damaged if the rotor is forced sideways or dropped onto the shaft.

- **2** Center the rotor over the drive shaft and carefully lower it straight down.
- **3** Secure the rotor to the drive shaft with the tie-down screw.
  - Tighten the screw by turning it to the right (clockwise) with the T-handle rotor wrench (365636 for both rotors).
- **4** Refer to the instrument instruction manual for centrifuge operation.
  - See Run Speeds, page 7, for information about speed limitations.

## **Removal and Sample Recovery**



If disassembly reveals evidence of leakage, you should assume that some fluid escaped the rotor. Apply appropriate decontamination procedures to the centrifuge and accessories.

- 1 Using the T-handle rotor wrench, turn the tie-down screw to the left (counterclockwise) to release the rotor.
- **2** Remove the rotor by lifting it straight up and off the drive shaft.

**3** Remove the rotor lid and remove tubes.

#### **Tubes**

The F301.5 and FX301.5 rotors each hold up to thirty 1.5 to 1.8-mL reaction-vials, listed in Table 1. The rotors also hold 250 to 400- $\mu$ L vials in adapters, commercially available 500 to 750- $\mu$ L tubes in adapters, as well as commercially available 2.0 and 2.2-mL tubes.

Refer to Chemical Resistances for information on the chemical compatibilities of labware materials.

Table 1 Available Tubes for the F301.5 and FX301.5 Rotors

Description	Dimensions	Volume	Part Number	Adapter	Max Speed (rpm)
polyethylene tube with attached cap	11 × 45 mm	1.8 mL	340196 (pkg/500)	a	10,000
polyallomer tube with attached cap (natural)	11 × 40 mm	1.5 mL	357448 (pkg/500)	a	14,000
polypropylene tube with separate cap	11 × 40 mm	1.5 mL	343169 (pkg/500)	a	14,000
polyethylene tube, plain	7 × 40 mm	400 μL	314326 (pkg/1000)	361247 (pkg/24)	11,500
polyethylene tube, Heparin- Lithium Fluoride coated	7 × 40 mm	400 μL	652824 (pkg/1000)	361247 (pkg/24)	11,500
polyethylene tube, Heparin- Lithium coated	7 × 40 mm	400 μL	652825 (pkg/1000)	361247 (pkg/24)	11,500
polyethylene tube, Heparin- Lithium Fluoride coated	5 × 45 mm	250 μL	652821 (pkg/1000)	361247 (pkg/24)	11,500
polyethylene tube, Heparin- Lithium coated	5 × 45 mm	250 μL	652822 (pkg/1000)	361247 (pkg/24)	11,500
polyethylene tube, plain	5 × 45 mm	250 μL	652823 (pkg/1000)	361247 (pkg/24)	11,500
500 to 750-μL tubes		500 μL/ 750 μL	commercially available	364690 (pkg/24)	11,500

a. No adapter is required with this tube; however, adapter 364701 (pkg/12) can be used.

1-6 MMR-TB-002DB



#### **Temperature Limits**

- Plastic containers have been centrifuge tested for use at temperatures between 2 and 25°C. For centrifugation at other temperatures, pretest containers under anticipated run conditions.
- If plastic containers are frozen before use, make sure that they are thawed to at least 2°C before centrifugation.

### **Run Speeds**

The centrifugal force at a given radius in a rotor is a function of speed. Comparisons of forces between different rotors are made by comparing the rotors' relative centrifugal fields (RCF). When rotational speed is adjusted so that identical samples are subjected to the same RCF in two different rotors, the samples are subjected to the same force.

The maximum run speed listed in the rotor specifications is for operation when all conditions are within the standard specifications. Speeds must be reduced under the following circumstances:

1. If nonprecipitating solutions more dense than 1.2 g/mL are centrifuged, the maximum allowable run speed must be reduced according to the following equation:

reduced maximum speed = (14,000 rpm\*) 
$$\sqrt{\frac{1.2 \text{ g/mL}}{\rho}}$$

\*Use 13,200 for nonrefrigerated Allegra X-22 centrifuges

where  $\rho$  is the density of the tube contents. This speed reduction will protect the rotor from excessive stresses due to the added tube load. Note, however, that the use of this formula may still produce maximum speed figures that are higher than the limitations imposed by the use of certain tubes or adapters. In such cases, use the lower of the two figures.


- **2.** Further speed limits must be imposed when self-forming-gradient salts are centrifuged, as the equation does not predict concentration limits/speeds that are required to avoid precipitation of salt crystals.
- **3.** Speeds may need to be reduced to maintain lower set temperatures (i.e. minimum temperature achievable at full rated speed in the Allegra X-30R is 8°C; to achieve a lower temperature the speed will need to be reduced).

MMR-TB-002DB 1-7

#### **Care and Maintenance**

#### **Maintenance**

Do not use sharp tools on the rotor, as they can scratch the anodized surface. Corrosion begins in scratches and may open fissures in the rotor with continued use.

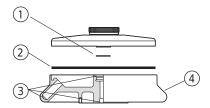


- 1. Lid O-ring (368896)
- 3. Check for Corrosion
- 2. O-ring (368991)
- 4. Rotor Body
- Periodically (at least monthly) inspect the rotor, especially inside cavities, for rough spots, cracks, pitting, white powder deposits (frequently aluminum oxide), or heavy discoloration.
  - If any of these signs are evident, do not run the rotor.
  - Contact your Beckman Coulter representative for information about the Field Rotor Inspection Program and the rotor repair center.
- **2** Before using the tie-down screw, check it for damage such as distortion, splitting, or stripped threads.
  - Replace it if it is damaged.
- **3** Regularly apply a thin, even coat of Spinkote lubricant (306812) to the rotor drive hole to prevent rotor sticking.

Refer to *Chemical Resistances* for the chemical compatibilities of rotor and accessory materials. Your Beckman Coulter representative provides contact with the Field Rotor Inspection Program and the rotor repair center.

1-8 MMR-TB-002DB

#### Cleaning

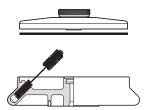

Wash the rotor and rotor components immediately if salts or other corrosive materials are used or if spillage has occurred. Do not allow corrosive materials to dry on the rotor.

Under normal use, wash the rotor frequently (at least weekly) to prevent buildup of residues. If the rotor is left in the centrifuge for long periods of time, remove it at least once a month for cleaning and lubrication.



- Remove the O-ring and wash the rotor, lid, and O-ring in a mild detergent, such as Beckman Solution 555 (339555), that won't damage the rotor.
  - The Rotor Cleaning Kit (339558) contains two plastic-coated brushes and two quarts of Solution 555 for use with rotors and accessories.
  - Dilute the detergent 10 to 1 with water.

**NOTE** Do not wash the rotor components or accessories in a dishwasher. Do not soak components in detergent solution for long periods of time, such as overnight.

- 2 Thoroughly rinse the cleaned rotor and components with distilled water.
- **3** Air-dry the rotor and lid upside down.
  - Do not use acetone to dry the rotor.
- 4 Lightly but evenly lubricate the rotor drive-hole threads with Spinkote (306812).
- 5 Apply a thin, even coat of silicone vacuum grease (335148) to the O-ring, then replace it in the groove in the outer rim of the rotor.
- **6** Clean metal threads every 6 months, or as necessary.
  - Use a brush and concentrated Solution 555.
  - Rinse and dry thoroughly, then lubricate lightly but evenly with Spinkote to coat all threads.

#### **Decontamination**





If aluminum rotor components become contaminated with radioactive material, decontaminate it using a solution that will not damage the anodized surfaces. Beckman Coulter has tested a number of solutions and found two that do not harm anodized aluminum: RadCon Surface Spray or IsoClean Solution (for soaking),\* and Radiacwash.†

**NOTE** IsoClean can cause fading of colored anodized surfaces. Use it only when necessary and remove it promptly from surfaces.

While Beckman Coulter has tested these methods and found that they do not damage the rotor or components, no guarantee of decontamination is expressed or implied. Consult your laboratory safety officer regarding the proper decontamination methods to use.

If the rotor or accessories are contaminated with toxic or pathogenic solutions, follow appropriate decontamination procedures. Check *Chemical Resistances* to be sure the decontamination method will not damage any part of the rotor.

#### Sterilization and Disinfection



- The rotor and all rotor components can be autoclaved at 121°C for up to one hour. Remove the lid from the rotor and place the rotor, lid, and O-ring in the autoclave upside down.
- Ethanol (70%)<sup>‡</sup> or hydrogen peroxide (6%) may be used on all rotor components, including those made of plastic. Bleach (sodium hypochlorite) may be used, but may cause discoloration of anodized surfaces. Use the minimum immersion time for each solution, per laboratory standards.

While Beckman Coulter has tested these methods and found that they do not damage the rotor or components, no guarantee of sterility or disinfection is expressed or implied. When sterilization or disinfection is a concern, consult your laboratory safety officer regarding proper methods to use.

Refer to publication IN-192, included in each box of tubes or bottles, for tube sterilization and disinfection procedures.

## **Storage**

When the rotor is not in use, store it in a dry environment (not in the centrifuge) with the rotor lid removed to allow air circulation so moisture will not collect in the tube cavities.

1-10 MMR-TB-002DB

<sup>\*</sup> In U.S., contact Nuclear Associates (New York); in Eastern Europe and Commonwealth States, contact Victoreen GmbH (Munich); in South Pacific, contact Gammasonics Pty. Ltd. (Australia); in Japan, contact Toyo Medic Co. Ltd. (Tokyo).

<sup>†</sup> In U.S., contact Biodex Medical Systems (Shirley, New York); internationally, contact the U.S. office to find the dealer closest to you.

<sup>‡</sup> Flammability hazard. Do not use in or near operating ultracentrifuges.

### **Returning a Rotor**

Before returning a rotor or accessory for any reason, prior permission must be obtained from Beckman Coulter, Inc. This form may be obtained from your local Beckman Coulter sales office. The form, entitled *Returned Material Authorization* (RMA) for United States returns or *Returned Goods Authorization* (RGA) for international returns, should contain the following information:

- rotor type and serial number,
- history of use (approximate frequency of use),
- reason for the return,
- original purchase order number, billing number, and shipping number, if possible,
- name and email address of the person to be notified upon receipt of the rotor or accessory at the factory,
- name and email address of the person to be notified about repair costs, etc.

To protect our personnel, it is the customer's responsibility to ensure that all parts are free from pathogens and/or radioactivity. Sterilization and decontamination must be done before returning the parts. Smaller items (such as tubes, bottles, etc.) should be enclosed in a sealed plastic bag.

All parts must be accompanied by a note, plainly visible on the outside of the box or bag, stating that they are safe to handle and that they are not contaminated with pathogens or radioactivity. **Failure to attach** this notification will result in return or disposal of the items without review of the reported problem.

Use the address label printed on the RGA/RMA form when mailing the rotor and/or accessories.

Customers located outside the United States should contact their local Beckman Coulter office.

### **Supply List**

**NOTE** Publications referenced in this manual can be obtained at www.beckman.com or by calling Beckman Coulter at 1-800-742-2345 in the United States, or by contacting your local Beckman Coulter office.

Call Beckman Coulter Sales (1-800-742-2345 in the United States) for detailed information on ordering parts and supplies. For your convenience, a partial list is given below.

MMR-TB-002DB 1-11

## **Replacement Rotor Parts**

F301.5 rotor assembly	368894
F301.5 lid assembly	368895
Tie-down screw, F301.5	365806
FX301.5 rotor assembly	392274
FX301.5 lid assembly	392275
Tie-down screw, FX301.5	361367
Lid O-ring	368896
Rotor O-ring	368991
T-handle rotor wrench	365636

**NOTE** The same lid O-ring, rotor O-ring, and T-handle wrench are used in both the F301.5 and FX301.5 rotors.

#### Other

Tubes and accessories	see Table 1
Spinkote lubricant (2 oz)	306812
Silicone vacuum grease (1 oz)	335148
Rotor Cleaning Kit	339558
Beckman Solution 555 (1 qt)	339555

1-12 MMR-TB-002DB

## Beckman Coulter, Inc. Benchtop Rotor Warranty

Subject to the conditions specified below and the warranty clause of the Beckman Coulter, Inc., terms and conditions in effect at the time of sale, Beckman Coulter agrees to correct either by repair or, at its election, by replacement, any defects of material or workmanship which develop within seven (7) years after delivery of a benchtop centrifuge rotor to the original buyer by Beckman Coulter or by an authorized representative, provided that investigation and factory inspection by Beckman Coulter discloses that such defect developed under normal and proper use. Should a Beckman Coulter centrifuge be damaged due to a failure of a rotor covered by this warranty, Beckman Coulter will supply free of charge all centrifuge parts required for repair.

#### **Conditions**

Except as otherwise specifically provided herein, this warranty covers the rotor only and Beckman Coulter shall not be liable for damage to accessories or ancillary supplies including but not limited to (i) tubes, (ii) tube caps, (iii) tube adapters, or (iv) tube contents.

This warranty is void if the rotor has been subjected to customer misuse such as operation or maintenance contrary to the instructions in the Beckman Coulter rotor or centrifuge manual.

This warranty is void if the rotor is operated with a rotor drive unit or in a centrifuge unmatched to the rotor characteristics or operated in a Beckman Coulter centrifuge that has been improperly disassembled, repaired, or modified.

Thermoplastic rotors or components used in some benchtop centrifuges are warranted for one (1) year from date of purchase.


#### Disclaimer

IT IS EXPRESSLY AGREED THAT THE ABOVE WARRANTY SHALL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND OF THE WARRANTY OF MERCHANTABILITY AND THAT NEITHER BECKMAN COULTER, INC. NOR ITS SUPPLIERS SHALL HAVE ANY LIABILITY FOR SPECIAL OR CONSEQUENTIAL DAMAGES OF ANY KIND WHATSOEVER ARISING OUT OF THE MANUFACTURE, USE, SALE, HANDLING, REPAIR, MAINTENANCE, OR REPLACEMENT OF THE PRODUCT.

MMR-TB-002DB Warranty-1

**Beckman Coulter, Inc. Benchtop Rotor Warranty** 

Warranty-2 MMR-TB-002DB

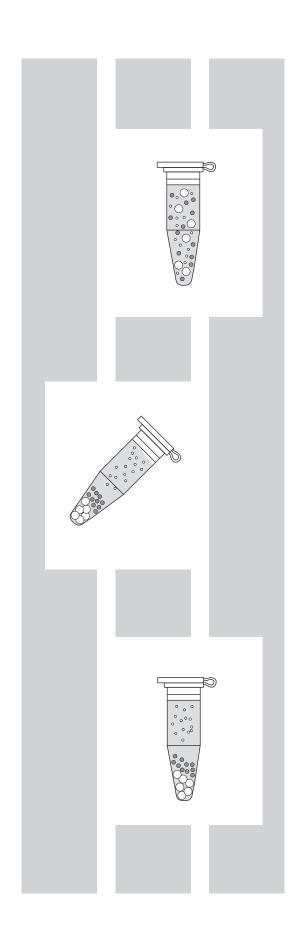

## Related Documents

#### Microfuge 22R (MMR-IM-5)

- Description
- Installation
- Operation
- Troubleshooting
- Care and Maintenance

#### Allegra X-22 Series (GS22-IM-6)

- Description
- Installation
- Operation
- Troubleshooting
- Care and Maintenance

#### Allegra X-30 Series (B01145)

- Safety
- Introduction
- Description
- Installation
- Operation
- Troubleshooting
- Care and Maintenance

Available in hard copy or electronic pdf by request.

#### **Additional References**

- Chemical Resistances for Beckman Coulter Centrifugation Products (IN-175)
- Use and Care of Tubes and Bottles (IN-192)

Available in hard copy or electronic pdf by request.

Available at www.beckman.com

www.beckman.com

