

JA-18 Fixed Angle Rotor

For Use in Beckman Coulter J2, J6, and Avanti J Series Centrifuges



PN J-TB-035AT November 2018





JA-18 Fixed Angle Rotor

PN J-TB-035AT (November 2018)

© 2018 Beckman Coulter, Inc. All rights reserved.

Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks mentioned herein are trademarks or registered trademarks of Beckman Coulter, Inc. in the United States and other countries.

All other trademarks, service marks, products, or services are trademarks or registered trademarks of their respective holders.

Contact Us

If you have any questions, contact our Customer Support Center.

- Worldwide, find us via our website at www.beckman.com/support/technical
- In the USA and Canada, call us at 1-800-369-0333.
- Outside of the USA and Canada, contact your local Beckman Coulter Representative.

Find us on the World Wide Web at: www.beckman.com

EC REP

Beckman Coulter Eurocenter S.A. 22, rue Juste-Olivier Case Postale 1044 CH - 1260 Nyon 1, Switzerland Tel: +41 (0) 22 365 36 11

Glossary of Symbols is available at beckman.com/techdocs (PN C24689).

Original Instructions

Revision History

For updates, go to www.beckman.com/techdocs and download the most recent manual or system help for your instrument.

Issue AT, 11/18 - Changes or additions were made in the following:

- Certified Free Tubes
- Sterile Tubes
- Table 1, Available Tubes and Bottles for the JA-18 Fixed Angle Rotor
- Sterilization and Disinfection

Changes that are part of the most recent revision are indicated in text by a bar in the left margin of the amended page.

PN J-TB-035AT iii

IV PN J-TB-035AT

Safety Notice

Read all product manuals and consult with Beckman Coulter-trained personnel before attempting to use this equipment. Do not attempt to perform any procedure before carefully reading all instructions. Always follow product labeling and manufacturer's recommendations. If in doubt as to how to proceed in any situation, contact your Beckman Coulter Representative.

Alerts for Warning, Caution, Important, and Note



WARNING indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury.



CAUTION indicates a potentially hazardous situation, which, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices.

IMPORTANT IMPORTANT is used for comments that add value to the step or procedure being performed. Following the advice in the Important adds benefit to the performance of a piece of equipment or to a process.

NOTE NOTE is used to call attention to notable information that should be followed during installation, use, or servicing of this equipment.

PN J-TB-035AT

Safety Information for the JA-18 Rotor

This safety notice summarizes information basic to the safe use of the rotor described in this manual. The international symbol displayed in *Alerts for Warning, Caution, Important, and Note* is a reminder to the user that all safety instructions should be read and understood before operation or maintenance of this equipment is attempted. When you see the symbol on other pages throughout this publication, pay special attention to the specific safety information presented. Observance of safety precautions will also help to avoid actions that could damage or adversely affect the performance of the rotor. This rotor was developed, manufactured, and tested for safety and reliability as part of a Beckman Coulter centrifuge/rotor system. Its safety or reliability cannot be assured if used in a centrifuge not of Beckman Coulter's manufacture or in a Beckman Coulter centrifuge that has been modified without Beckman Coulter's approval.

! WARNING

Handle body fluids with care because they can transmit disease. No known test offers complete assurance that such fluids are free of micro-organisms. Some of the most virulent — Hepatitis (B and C) viruses, HIV (I -V), atypical mycobacteria, and certain systemic fungi — further emphasize the need for aerosol protection. Handle other infectious samples according to good laboratory procedures and methods to prevent spread of disease. Because spills may generate aerosols, observe proper safety precautions for aerosol containment. Do not run toxic, pathogenic, or radioactive materials in this rotor without taking appropriate safety precautions. Biosafe containment should be used when Risk Group II materials (as identified in the World Health Organization Laboratory Biosafety Manual) are handled; materials of a higher group require more than one level of protection.

MARNING

The rotor and accessories are not designed for use with materials capable of developing flammable or explosive vapors. Do not centrifuge such materials in nor handle or store them near the centrifuge.

MARNING

Although rotor components and accessories made by other manufacturers may fit in the JA-18 rotor, their safety in this rotor cannot be ascertained by Beckman Coulter. Use of other manufacturers' components or accessories in the JA-18 rotor may void the rotor warranty and should be prohibited by your laboratory safety officer. Only the components and accessories listed in this publication should be used in this rotor.

vi PN J-TB-035AT

⚠ WARNING

Do not run an empty rotor. Place filled tubes in at least two opposing cavities. Make sure that filled containers are loaded symmetrically into the rotor and that opposing tubes are filled to the same level with liquid of the same density.

MARNING

If disassembly reveals evidence of leakage, and pathogenic or radioactive materials are involved, you should assume that some fluid escaped the rotor. Apply appropriate decontamination procedures to the centrifuge and accessories.

MARNING

Never exceed the maximum rated speed of the rotor and labware in use. Refer to the section on Run Speeds.

A CAUTION

Do not use sharp tools on the rotor that could cause scratches in the rotor surface. Corrosion begins in scratches and may open fissures in the rotor with continued use.

PN J-TB-035AT vii

Safety Notice Safety Information for the JA-18 Rotor

PN J-TB-035AT viii

Contents

```
Revision History, iii
Safety Notice, v
Alerts for Warning, Caution, Important, and Note, v
Safety Information for the JA-18 Rotor, vi
JA-18 Fixed Angle Rotor, 1
Specifications, 1
Description, 2
Preparation and Use, 3
         Prerun Safety Checks, 3
         Rotor Preparation, 3
         Tubes and Bottles, 4
            Temperature Limits, 4
             Certified Free Tubes, 4
            Sterile Tubes, 5
            Thickwall Tubes, 5
            Polycarbonate and Polypropylene Bottles, 5
             Available Tubes and Bottles for the JA-18 Fixed Angle Rotor, 5
         Operation, 8
         Installing the Rotor, 8
         Removal and Sample Recovery, 10
         Temperature, 12
             Avanti J Series Centrifuges, 12
             Microprocessor-Controlled J2 and J6 Series Centrifuges
                (Models J2-MI, J2-MC, J6-MI, and J6-MC), 12
             Analog J2 Series Centrifuges, 14
         Run Times, 15
         Run Speeds, 16
Care and Maintenance, 17
         Maintenance, 17
         Cleaning, 18
         Decontamination, 19
         Sterilization and Disinfection, 19
         Storage, 20
Returning a Rotor, 20
Supply List, 21
         Replacement Rotor Parts, 21
         Other, 21
```

Beckman Coulter, Inc. J Series Rotor Warranty Related Documents

Tables


1	Available Tubes and Bottles for the JA-18 Fixed Angle Rotor, 5
2	Temperature Compensation Settings — Models J2-MI and J2-MC Centrifuges, 13
3	${\it Temperature Compensation Settings-J6 Series Centrifuges, 13}$
4	Temperature Compensation Settings — Models J2-HS and J2-21 Centrifuges, 14
5	Temperature Compensation Settings — Model J2-HC Centrifuge, 14
6	Relative Centrifugal Fields for the JA-18 Rotor, 16

Illustrations

1	Arranging Tubes or Bottles in the Rotor, 3
2	Rotor Drive Pin Location and Orientation, 9
3	Centrifuge Drive Spindle Hub Configurations, 9
4	Using the Rotor Removal Tool, 11

JA-18 Fixed Angle Rotor

Specifications

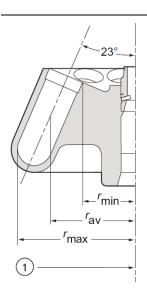

1. Axis of Rotation

Maximum speed	18,000 rpm
Critical speed range ^a	600 to 800 rpm
Maximum solution density	1.2 g/mL
Maximum allowable imbalance of opposing loads	2 grams
Relative Centrifugal Field ^b at maximum speed	
At <i>r_{max}</i> (132 mm)	47,900× <i>g</i>
At r _{av} (98 mm)	35,600× <i>g</i>
At <i>r_{min}</i> (64 mm)	23,200×g
k factor at maximum speed	566
Number of tube cavities	10
Available tubes and bottles	see Table 1
Nominal dimensions of largest tube	38×102 mm
Nominal tube capacity	100 mL
Nominal rotor capacity	1 L
Approximate acceleration time to maximum speed	
(rotor fully loaded)	6.5 min
Approximate deceleration time from maximum speed	
(rotor fully loaded, max. brake)	5 min
Weight of fully loaded rotor	12 kg (26 lb)
Rotor material	aluminum
Conditions requiring speed reduction	see Run Speeds
Rotor entry code	


- a. The critical speed range is the range of speeds over which the rotof shifts so as to rotate about its center mass. Passing through the critical speed range is characterized by some vibration.
- b. Relative Centrifugal Field (RCF) is the ratio of the centrifugal acceleration at a specified radius and speed $(r\omega^2)$ to the standard acceleration of gravity (g) according to the following formula:

$$RCF = \frac{r\omega^2}{g}$$

where r is the radius in millimeters, ω is the angular velocity in radians per second (2 π RPM /60), and g is the standard acceleration of gravity (9807 mm/s²). After substitution:

$$RCF = 1.12r \left(\frac{RPM}{1000}\right)^2$$

Description

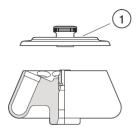

1. Lid (PN 364845)

This rotor has been manufactured in a registered ISO 9001 or 13485 facility for use with the specified Beckman Coulter centrifuges.

The JA-18 fixed angle rotor, rated for 18,000 rpm, is designed to centrifuge up to ten 100-mL containers at a 23-degree angle to the axis of rotation. Used in Beckman Coulter J2, J6, and Avanti J series centrifuges, the rotor develops centrifugal forces that can efficiently pellet cells, cellular organelles, and membrane fragments.

The rotor body and lid are made of aluminum and are anodized for corrosion resistance — the rotor is black and the lid is blue.

Pins in the rotor drive hole mate with grooves in the centrifuge drive hub (older model centrifuges), or with drive hub teeth (new model centrifuges), to prevent the rotor from slipping during acceleration and deceleration.

Starting in 2002, JA-18 rotors are manufactured with magnets embedded in the rotor base for rotor identification in the Avanti J-E centrifuge. Older JA-18 rotors without magnets can be used in the Avanti J-E, but the rotor speed will be limited to 14,000 rpm. (Other Avanti J series and J2 series centrifuges use alternate methods of rotor identification that enable the rotor to run up to 18,000 rpm.)

The rotor has been tested 1 to demonstrate containment of microbiological aerosols under normal operating conditions of the associated Beckman Coulter centrifuge, when used and maintained as instructed.

The JA-18 rotor is warranted for 7 years (see the Beckman Coulter, Inc. J Series Rotor Warranty).

1-2 PN J-TB-035AT

¹ Validation of microbiological containment was done at an independent third-party testing facility (CAMR, Porton Down, UK, or USAMRIID, Ft. Detrick, MD, U.S.A.). Improper use or maintenance may affect seal integrity and thus containment.

Preparation and Use

Specific information about the JA-18 rotor is given here. Use the Rotors and Tubes for Beckman Coulter J2, J6, and Avanti J Series Centrifuges (PN JR-IM-10) along with this rotor manual for complete rotor and accessory operation.

NOTE Although rotor components and accessories made by other manufacturers may fit in the JA-18 rotor, their safety in this rotor cannot be ascertained by Beckman Coulter. Use of other manufacturers' components or accessories in this rotor may void the rotor warranty and should be prohibited by your laboratory safety officer. Only the components and accessories listed in this publication should be used in this rotor.

Prerun Safety Checks

IMPORTANT Read the *Safety Notice* before using the rotor.


- 1 Make sure that the rotor and lid are clean and show no signs of corrosion or cracking
- **2** Check the chemical compatibilities of all materials used (refer to *Appendix A* in *Rotors and Tubes for Beckman Coulter J2, J6, and Avanti J Series Centrifuges* (PN JR-IM-10).
- **3** Verify that the tubes and bottles being used are listed in Table 1.

Rotor Preparation

For runs at other than room temperature, refrigerate or warm the rotor beforehand for fast temperature equilibration.

Load the filled containers symmetrically into the rotor. (Refer to *Tubes and Bottles* on page 1-4 for information about containers.) If fewer than ten tubes are being run, they must be arranged symmetrically in the rotor (see Figure 1). Opposing tubes must be filled to the same level with liquid of the same density (within 2 grams).

Figure 1 Arranging Tubes or Bottles in the Rotor

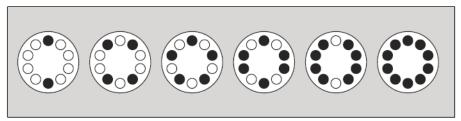

Two, four, five, six, eight, or ten tubes or bottles can be run, if they are arranged symmetrically in the rotor, as shown.

PN J-TB-035AT 1-3

Tubes and Bottles

The JA-18 rotor uses the tubes and bottles listed in Table 1. Use only those items listed, and observe the maximum speed limits and fill volumes shown. (Maximum fill volume is the maximum amount that can be centrifuged in the container listed.) Refer to Appendix A in Rotors and Tubes for Beckman Coulter J2, J6, and Avanti J Series Centrifuges (PN JR-IM-10) for chemical compatibilities of tube, bottle, and accessory materials.

Some conical tubes or bottles from other manufacturers may be used in this rotor with adapters. Because Beckman Coulter cannot test all manufacturers' tubes or bottles with respect to their ability to withstand the *g*-forces generated by this rotor, it is highly recommended that you pretest any other manufacturer's tubes in these adapters using water instead of valuable samples.



Some tube/adapter combinations (15-mL and 50-mL conical tubes with adapters PN 356592 and PN 356593; see Table 1) are too tall to allow the lid to seat properly on the rotor, so the lid cannot be used when these tubes are run. In Avanti J series centrifuges, the lid must be used during each run, so these tube/adapter combinations cannot be run in an Avanti J. In J2 and J6 series centrifuges, the rotor can be run without the lid; however, the rotor may not reach top speed and temperature control may also be impaired.



Do not run 50-mL tubes with snap-on caps in this rotor. The g-forces generated during operation may cause the snap-on caps to come off, which could dislodge the rotor lid.

Temperature Limits



- Plastic containers have been centrifuge tested for use at temperatures between 2 and 25°C. For centrifugation at other temperatures, pretest tubes under anticipated run conditions.
- If plastic containers are frozen before use, make sure that they are thawed to at least 2°C prior to centrifugation.

Certified Free Tubes



Certified free tubes are lot traceable to testing that confirms the absence of endotoxin, DNase, RNase, and human & mouse DNA below a detectable limit.


1-4 PN J-TB-035AT

Sterile Tubes



Sterile tubes are sterilized via ethylene oxide in compliance with ISO 11135:2014. Cartons include several peel packages, each containing a typical run quantity of tubes. Packaging meets requirements of ISO11607:2006.

Thickwall Tubes

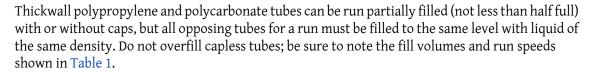

Polycarbonate and Polypropylene Bottles

Capped polycarbonate and polypropylene bottles may be run completely filled, or partially filled (not less than half full). All opposing containers for a run must be filled to the same level with liquid of the same density. Be sure to note the fill volumes and run speeds shown in Table 1.

Available Tubes and Bottles for the JA-18 Fixed Angle Rotor

Table 1 Available Tubes and Bottles for the JA-18 Fixed Angle Rotor^a

	Tube	Required A				
Dimensions and Volume	Description	Part Number	Max. Fill Volume (mL)	Description	Part Number	Max Speed ^b / RCF/ k factor
	Certified Free & Sterile Polypropylene Quick Seal	C14305 Carton of 24 (4 packs of 6)		none	_	18,000 rpm 47,900 × <i>g</i> 566
38 × 102 mm 100 mL	Certified Free Polypropylene Quick Seal	C14289 (pkg/25)	100			
	Standard Polypropylene Quick Seal	345776 (pkg/25)				





 Table 1
 Available Tubes and Bottles for the JA-18 Fixed Angle Rotor^a (Continued)

	Tube			Required A		
Dimensions and Volume	Description	Part Number	Max. Fill Volume (mL)	Description	Part Number	Max Speed ^b / RCF/ <i>k</i> factor
38 × 102 mm 100 mL	Ultra-Clear Quick-Seal tube	345778	100	none	_	18,000 rpm 47,900 × <i>g</i> 566
38 × 102 mm 100 mL	polypropylene bottle w/cap assembly	355624 ^c	94	none	_	18,000 rpm 47,900 × <i>g</i> 566
38 × 102 mm 100 mL	polycarbonate bottle w/cap assembly	355620	70	none	_	18,000 rpm 47,900 × <i>g</i> 566
29 × 104 mm 50 mL	polycarbonate open-top tube ^d	363647	50	adapter	347539 ^e	18,000 rpm 44,000 × g —
29 × 104 mm 50 mL	polycarbonate open-top tube ^d	357007	50	adapter	347539 ^e	18,000 rpm 44,000 × <i>g</i> —
29×104 mm 50 mL	polycarbonate bottle w/screw cap	357002	34	adapter	347539 ^e	18,000 rpm 44,000 × <i>g</i> —
29×104 mm 50 mL	polypropylene bottle w/screw cap	357003	34	adapter	347539 ^e	18,000 rpm 44,000 × <i>g</i> —
11 × 38 mm 1.8 mL	polyethylene microcentrifuge tube	340196	1.8	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
11 × 39 mm 1.5 mL	polypropylene microcentrifuge tube (natural)	357448	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
11 × 39 mm 1.5 mL	polypropylene microcentrifuge tube (natural)	356090	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
11 × 39 mm 1.5 mL	polypropylene microcentrifuge tube (blue)	356091	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
11 × 39 mm 1.5 mL	polypropylene microcentrifuge tube (green)	356092	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —

1-6 PN J-TB-035AT

Table 1 Available Tubes and Bottles for the JA-18 Fixed Angle Rotor^a (Continued)

	Tube		Required A			
Dimensions and Volume	Description	Part Number	Max. Fill Volume (mL)	Description	Part Number	Max Speed ^b / RCF/ <i>k</i> factor
11 × 39 mm 1.5 mL	polypropylene microfuge tube (yellow)	356093	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
11 × 39 mm 1.5 mL	polypropylene microcentrifuge tube (orange)	356094	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × g —
11 × 38 mm 1.5 mL	polypropylene microcentrifuge tube (natural) (separate cap)	343169	1.5	adapter	344497 ^h inside 347539 ^e	18,000 rpm 35,000 × <i>g</i> —
30 × 115 mm 50 mL	conical tube	f	37	adapter	356963 ^g h	14,000 rpm 29,000 × <i>g</i> —
17 × 120 mm 15 mL	conical tube	f	15	adapter	356963 ^g h	15,000 rpm ⁱ 33,300 × <i>g</i> —
13 × 100 mm 10 mL	round-bottom tube	f	10	adapter	356975 ^h j	14,000 rpm ⁱ 29,000 × <i>g</i> —
12 × 75 mm 3 mL	round-bottom tube	f	3	adapter	356969 ^h j	14,000 rpm ⁱ 29,000 × <i>g</i> —

- a. Use only the items listed here and observe fill volumes and maximum run speeds.
- b. Maximum speeds listed are guidelines only. These speeds have been achieved in reliability tests at Beckman Coulter, but, because of manufacturing variances, no guarantee of performance or fit is expressed or implied
- c. This part number is for the complete bottle assembly; to order separate components, part numbers are: bottle only, PN 355626; cap, PN 355625; cap O-ring, PN 889633
- d. A cap is available for these tubes, but it cannot be used in the JA-18 rotor. For a cap to use when storing samples, order cap PN 356262.
- e. Polypropylene.
- f. Commercially available. Observe manufacturer's speed and fill volume guidelines.
- g. Do not use the rotor lid when using these tube/adapter combinations, as the lid will not seat properly. Do not use these tube/adapter combinations in this rotor in an Avanti J series centrifuge.
- h. Delrin.
- i. Recommended speed at temperature of 20°C or less; at higher temperatures, reduce speed to 11,000 rpm.
- j. Each adapter holds three tubes.

PN J-TB-035AT 1-7

Operation

- 1 For low-temperature runs, precool the rotor in the centrifuge or in a refrigerator before use especially before short runs to ensure that the rotor reaches the set temperature. A suggested precooling cycle is a minimum of 30 minutes at 2000 rpm at the required temperature.
- 2 If fluid containment is required, use capped tubes or bottles. It is strongly recommended that all containers carrying physiological fluids be capped to prevent leakage.
- **3** If you are using an Avanti J series centrifuge (except Avanti J-E), select the JA-18.
- 4 If you are using an Avanti J-E or a microprocessor-controlled J2 or J6 series centrifuge, enter rotor code 18. Refer to your centrifuge instruction manual for more information.

Installing the Rotor

1 Lubricate J2 and J6 series centrifuge drive hubs with a thin, even coat of Spinkote (PN 306812) to prevent the rotor from sticking.


NOTE Avanti J series centrifuges have Delrin rings on the drive hubs to prevent rotor sticking, and do not require lubrication.


TIP Prior to installing the rotor, consider the following:

- Two metal pins inside the rotor drive hole engage with teeth on the centrifuge drive spindle hub to prevent the rotor from slipping during acceleration and deceleration (see Figure 2). When the rotor is properly installed, the pins sit either in the drive hub grooves or on top of the hub surface (see Figure 3).
- The rotor pins are positioned parallel to the BECKMAN COULTER name engraved at the center
 of the rotor body. Knowing the pin orientation before you install the rotor will help to ensure that
 you position the rotor properly on the hub, minimizing the chance of hub damage.

1-8 PN J-TB-035AT


Figure 2 Rotor Drive Pin Location and Orientation

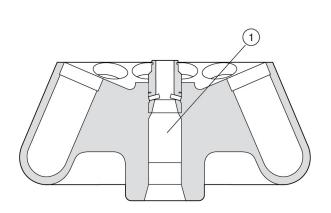

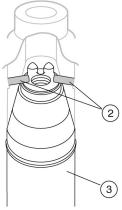

- 1. Rotor Drive Hole
- 2. Drive Pins (Angled pins shown. Pins can also be vertical or horizontal.
- 3. Avanti J Centrifuge Drive Spindle

Figure 3 Centrifuge Drive Spindle Hub Configurations

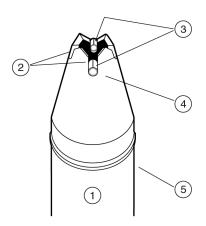

1. Older Model Centrifuges — Be sure the pins

in the rotor drive hole are seated in the


- (6)
- grooves of the drive spindle hub as shown. 2. Drive Hub Grooves (4)
- 3. Rotor Pins (2)
- 4. Drive Hub
- 5. Drive Spindle Assembly

6. Newer Model Centrifuges — Be sure the pins in the rotor drive hole are not sitting on top of the teeth on the drive spindle hub.


(9)

- 7. Teeth (2)
- 8. Drive Hub
- 9. Drive Spindle Assembly
- Turn the rotor until the BECKMAN COULTER name is either horizontal or perpendicular to the drive hub pins or teeth. Alternately, rotate the drive spindle to align the hub pins or teeth with the rotor drive pins.

PN J-TB-035AT 1-9

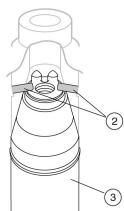

- **3** Carefully lower the rotor straight down onto the centrifuge drive spindle hub. Do not drop the rotor onto the hub.
- 4 Slowly turn the rotor, making sure that it is properly seated on the drive spindle hub. In rare cases, the rotor pins can rest on top of the hub pins or teeth and not seat on the hub surface. This can prevent the rotor from being secured to the drive spindle and may cause a rotor mishap during operation.
- Place the lid on the rotor, fitting the underside of the lid knob over the hex bolt. The lid knob acts as a wrench for tightening and loosening the tie-down mechanism.

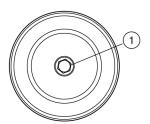

1. Underside of Lid Knob (use to tighten hex bolt)

6 Turn the lid knob to the right (clockwise) to attach the rotor to the hub.

Removal and Sample Recovery




If disassembly reveals evidence of leakage, you should assume that some fluid escaped the rotor. Apply appropriate decontamination procedures to the centrifuge and accessories.

1 Unscrew the rotor lid to release the rotor from the spindle hub.

1-10 PN J-TB-035AT

- Remove the rotor by lifting it straight up and off the drive spindle. If the rotor sticks to the drive spindle, a rotor removal tool (PN 346965) may be used (see Figure 4).
 - **a.** Remove the lid and use an adjustable wrench to unscrew the tie-down bolt assembly from the rotor. Set the bolt assembly aside.
 - **b.** Screw the larger end of the rotor removal tool into the threaded opening. As the removal tool tightens down, it pushes against the drive hub, forcing the rotor up and off of the hub.
 - **c.** Unscrew the removal tool.
 - **d.** Reassemble the tie-down bolt in the rotor and tighten it securely with the adjustable wrench.
 - **e.** Remove the rotor from the centrifuge.

Figure 4 Using the Rotor Removal Tool

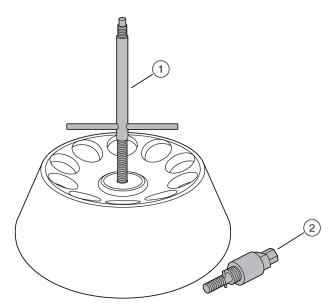

- 1. Rotor Removal Tool (PN 346965)
- 2. Tie-Down Bolt

PN J-TB-035AT 1-11

Temperature

To ensure that the JA-18 rotor reaches the required temperature during the run, follow the appropriate instructions below for the model of J centrifuge being used.

Avanti J Series Centrifuges

Enter the run temperature according to the instructions in your centrifuge instruction manual. No additional input is required.

Microprocessor-Controlled J2 and J6 Series Centrifuges (Models J2-MI, J2-MC, J6-MI, and J6-MC)

Enter rotor code 18; then follow the steps below.

- 1 Press the **[TEMP]** key on the centrifuge control panel, and then use the keypad to enter the sample temperature.
- **2** Find the compensation value in Table 2 or Table 3 that corresponds with the set temperature and run speed. The compensation values listed are approximates for all centrifuge models. If precise temperature control is required, determine the compensation settings empirically as follows:
 - **a.** Fill the rotor with water-filled tubes or bottles. Precool the rotor and water to the required temperature.
 - **b.** Place the filled, precooled rotor in the centrifuge and run the system for at least 1/2 hour.
 - **c.** Measure the water temperature. If it is lower than the required run temperature, enter positive compensation units during run setup following steps 3 and 4 below. If it is higher than the required run temperature, enter negative compensation units during run setup, also using steps 3 and 4 below.
 - **d.** Discard the water. Fill tubes with sample, load tubes into the rotor, and precool the system to the required temperature before starting the run.
- **3** Press [COMP ADJ]. the word "COMP" flashes below the TEMPERATURE display, and the display flashes.
- 4 Use the keypad to enter the desired compensation value. Press the [±.] key to enter a minus sign; pressing it again will remove the minus sign.
- **5** Check the temperature display. (If the entry is incorrect, press **[CE]** and reenter the digits.)

1-12 PN J-TB-035AT

6 When the entry is correct, press [ENTER/RECALL].

NOTE To clear a COMP ADJ entry, press [COMP ADJ], [0], and [ENTER/RECALL].

Table 2 Temperature Compensation Settings^a — Models J2-MI and J2-MC Centrifuges

Rotor Speed	Required Sample Temperature ^b (°C)							
(rpm)	-20°	-10°	2°	5°	10°	20°	40°	
18,000	N	N	N	N	N	-9	-7	
17,000	N	N	N	-9	-9	-8	-6	
16,000	N	N	–9	-9	-9	-8	-5	
12,000	N	-6	-8	-6	-4	-4	-3	
8,000	-5	-4	-3	-3	-2	-1	0	
5,000	-4	-3	-2	-2	-1	0	0	

a. Interpolate if intermediate values are required.

Table 3 Temperature Compensation Settings^a — J6 Series Centrifuges

Rotor Speed	Required Sample Temperature ^b (°C)					
(rpm)	2°	10°	20°	40°		
6,000	-3	-2	0	N		
4,000	-2	0	0	N		
2,000	-1	0	0	N		

a. Interpolate if intermediate values are required.

PN J-TB-035AT 1-13

b. An "N" indicates that the rotor cannot achieve the required temperature at this speed.

b. An "N" indicates that the rotor cannot achieve the required temperature at this speed.

Analog J2 Series Centrifuges

Enter the required run temperature and the appropriate temperature compensation units (see Table 4 or Table 5) on the centrifuge control panel as follows.

- 1 Turn the **SET** knob to the required sample temperature.
- **2** Find the compensation value that corresponds to the required temperature and run speed. The compensation values listed are approximate for all centrifuge models. If precise temperature control is required, follow the steps in *Microprocessor-Controlled J2 and J6 Series Centrifuges* (Models J2-MI, J2-MC, J6-MI, and J6-MC).

Table 4 Temperature Compensation Settings^a — Models J2-HS and J2-21 Centrifuges

Rotor Speed	Required Sample Temperature ^b (°C, green bar)							
(rpm)	-20°	-10°	2°	5°	10°	20°	40°	
18,000	N	N	N	N	N	-10	-7	
17,000	N	N	N	-10	-9	-8	-6	
16,000	N	N	-10	-10	-9	-7	-5	
12,000	N	-6	-6	-6	-5	-4	-3	
8,000	-5	-4	-4	-3	-2	-1	0	
5,000	-4	-3	-2	-2	-1	0	0	

a. Interpolate if intermediate values are required.

Table 5 Temperature Compensation Settings^a — Model J2-HC Centrifuge

Rotor Speed	Required Sample Temperature ^b (°C)							
(rpm)	-20°	-10°	2°	5°	10°	20°	40°	
18,000	N	N	N	N	N	-8	-4	
17,000	N	N	N	-10	-9	-8	-3	
16,000	N	N	-10	-9	-8	-7	-2	
12,000	N	-6	-6	-5	-4	-3	–1	
8,000	-3	-2	-2	-1	-1	0	N	
5,000	-1	-1	-1	-1	-0	+1	N	

a. Interpolate if intermediate values are required.

1-14 PN J-TB-035AT

b. An "N" indicates that the rotor cannot achieve the required temperature at this speed.

b. An "N" indicates that the rotor cannot achieve the required temperature at this speed.

Run Times

The k factor of the rotor is a measure of the rotor's pelleting efficiency. Beckman Coulter has calculated the k factors for all of its rotors, at maximum speed with full tubes, using the following formula:

$$k = \frac{\ln(r_{\text{max}}/r_{\text{min}})}{\omega^2} \times \frac{10^{13}}{3600}$$
 EQ 1

where ω is the angular velocity of the rotor in radians per second (ω = 0.105 × rpm), r_{max} is the maximum radius, and r_{min} is the minimum radius.

After substitution:

$$k = \frac{(2.533 \times 10^{11}) \ln(r_{\text{max}} / r_{\text{min}})}{\text{rpm}^2}$$

In the JA-18 rotor, the k factor is 566 for full 100-mL tubes at maximum speed. For containers that nominally hold less than 100 mL, the k factors will be different. Calculate the k factors for these containers using equation **EQ 2**.

Use the k factor in the following equation to estimate the run time t (in hours) required to pellet particles of known sedimentation coefficient s (in Svedberg units, S).

$$t = \frac{k}{s}$$
 EQ 3

Run times can be estimated for centrifugation at less than maximum speed by adjusting the k factor as follows:

$$k_{\text{adj}} = k_{\text{rated}} \left(\frac{18,000}{\text{actual run speed}} \right)^2$$

Run times can also be estimated from data established in prior experiments using a different rotor if the k factor of the previous rotor is known. For any two rotors, a and b,

$$\frac{t_{\rm a}}{t_{\rm b}} = \frac{k_{\rm a}}{k_{\rm b}}$$
 EQ 5

where the *k* factors have been adjusted for the actual run speed used.

NOTE For more information on *k* factors, see www.beckman.com/centrifuges/rotors/calculator.

PN J-TB-035AT 1-15

Run Speeds

The centrifugal force at a given radius in a rotor is a function of speed. Comparisons of forces between different rotors are made by comparing the rotors' relative centrifugal fields (RCF). When rotational speed is adjusted so that identical samples are subjected to the same rcf in two different rotors, the samples are subjected to the same force. The RCF at a number of rotor speeds is provided in Table 6.

Do not select rotational speeds higher than the maximums you have determined to be appropriate for your protocols (in no case above 18,000 rpm or the limit shown in Table 1). In addition, when centrifuging solutions more dense than 1.2 g/mL in this rotor, calculate the reduced maximum allowable rotor speed:

reduced maximum speed = (18,000 rpm)
$$\sqrt{\frac{1.2 \text{ g/mL}}{\text{density of tube contents}}}$$

Note, however, that the use of this formula may still produce maximum speed figures higher than the limitations imposed by the use of certain tubes or adapters (see Table 1). In such cases, use the lower of the two figures.


Table 6 Relative Centrifugal Fields for the JA-18 Rotor^a

Batan Guard	Relative Centrifugal Field (x g)						
Rotor Speed	At r _{max}	At r _{av}	At r _{min}				
(rpm)	(132 mm)	(98 mm)	(64 mm)				
18,000	47,900	35,600	23,200				
17,000	42,700	31,700	20,700				
16,000	37,800	28,100	18,400				
15,000	33,300	24,700	16,100				
14,000	29,000	21,500	14,000				
13,000	25,000	18,500	12,100				
12,000	21,300	15,800	10,300				
11,000	17,900	13,300	8,670				
10,000	14,800	11,000	7,170				
9,000	12,000	8,890	5,810				
8,000	9,460	7,020	4,590				
7,000	7,240	5,380	3,510				
6,000	5,320	3,950	2,580				
5,000	3,700	2,740	1,790				
4,000	2,370	1,760	1,150				
3,000	1,330	987	645				
2,000	591	439	286				
1,000	147	109	71				

a. Entires in this table are calculated from the formula: $RCF = 1.12 \ r \ (RPM/1000)^2$ and then rounded to three significant digits.


1-16 PN J-TB-035AT

Relative Centrifugal Fields, JA-18 Rotor

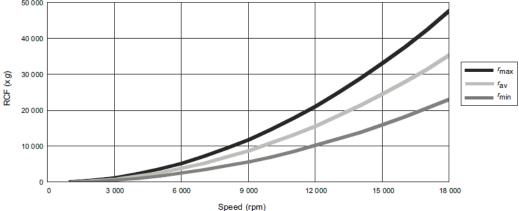

Care and Maintenance

Maintenance

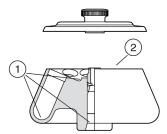

- 1. Check for Corrosion
- 2. Rotor Body

NOTE Do not use sharp tools on the rotor that could cause scratches in the rotor surface. Corrosion begins in scratches and may open fissures in the rotor with continued use.

- Periodically (at least monthly) inspect the rotor, especially inside cavities, for rough spots or
 pitting, white powder deposits (frequently aluminum oxide), or heavy discoloration. If any of
 these signs are evident, do not run the rotor. Contact us for information about the Field Rotor
 Inspection Program and the rotor repair center.
- Before using the rotor, inspect the rotor drive pins to ensure that they are not damaged. Damaged drive pins can prevent the rotor from seating properly on the centrifuge drive hub (Avanti J series centrifuges only). To inspect the drive pins, turn the rotor upside down and look into the drive hole in the center of the rotor. If the drive pins appear damaged, contact us.
- Regularly lubricate the metal threads in the rotor with a thin, even coat of Spinkote lubricant (PN 306812). Failure to keep these threads lubricated can result in damaged threads.
- Regularly apply Spinkote lubricant (PN 306812) to the centrifuge drive spindle to prevent the rotor from sticking (not required with Avanti J series centrifuges).

PN J-TB-035AT 1-17

Refer to Appendix A in Rotors and Tubes for Beckman Coulter J2, J6, and Avanti J Series Centrifuges (PN JR-IM-10) for chemical compatibilities of rotor and accessory materials. Your Beckman Coulter representative provides contact with the Field Rotor Inspection Program and the rotor repair center.

Cleaning


Wash the rotor and rotor components immediately if salts or other corrosive materials are used or if spillage has occurred. Do not allow corrosive materials to dry on the rotor.

Under normal use, wash the rotor frequently (at least weekly) to prevent buildup of residues.

1 Wash the rotor and lid in a mild detergent, such as Beckman Solution 555 (339555), that won't damage the rotor.

The Rotor Cleaning Kit (PN 339558) contains two special plastic-coated brushes and two quarts of Solution 555 for use with rotors and accessories.

Dilute the detergent 10 to 1 with water.

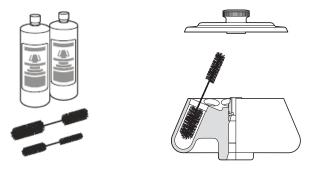

NOTE Do not wash rotor components in a dishwasher. Do not soak in detergent solution for long periods, such as overnight.

- **2** Thoroughly rinse the cleaned rotor and components with distilled water.
- **3** Air-dry the rotor and lid upside down. Do not use acetone to dry the rotor.

Clear metal threads every 6 months or as necessary. Use a brush and concentrated Solution 555. Rinse and dry thoroughly, then lubricate lightly but evenly with Spinkote to coat all threads.

NOTE The one-piece cap assembly for polypropylene bottle assembly PN 355624 contains an O-ring seal. If you remove the O-ring for cleaning or replacement, be careful not to damage the groove in the cap. The bottle cap and sealing surface of the O ring must be thoroughly dry before use.

1-18 PN J-TB-035AT

Decontamination





If the rotor (and/or accessories) becomes contaminated with radioactive material, it should be decontaminated using a solution that will not damage the anodized surfaces. Beckman Coulter has tested a number of solutions and found two that do not harm anodized aluminum: RadCon Surface Spray or IsoClean Solution (for soaking)¹, and Radiacwash².

NOTE IsoClean can cause fading of colored anodized surfaces. Use it only when necessary and remove it promptly from surfaces.

While Beckman Coulter has tested these methods and found that they do not damage components, no guarantee of decontamination is expressed or implied. Consult your laboratory safety officer regarding the proper decontamination methods to use.

If the rotor or other components are contaminated with toxic or pathogenic materials, follow appropriate decontamination procedures as outlined by your laboratory safety officer.

Sterilization and Disinfection



• The rotor and all rotor components can be autoclaved at 121°C for up to an hour. Remove the lid from the rotor and place the rotor and lid in the autoclave upside down.



Risk of injury or equipment damage. Ethanol is a flammability hazard. Do not use it in or near operating centrifuges.

• Ethanol (70%)³ may be used on all rotor components, including those made of plastic. Bleach (sodium hypochlorite) may be used, but may cause discoloration of anodized surfaces. Use the minimum immersion time for each solution, per laboratory standards.

While Beckman Coulter has tested these methods and found that they do not damage the rotor or components, no guarantee of sterility or disinfection is expressed or implied. When sterilization or disinfection is a concern, consult your laboratory safety officer regarding proper methods to use.

Where sterilization is critical in your application, consider using Beckman Coulter Certified Free & Sterilized Tubes. For tubes not available in the sterilized option, refer to publication IN-192, which is included in each box of tubes or bottles, for sterilization and disinfection procedures.

PN J-TB-035AT

¹ In U.S., contact Nuclear Associates (New York); in Eastern Europe and Commonwealth States, contact Victoreen GmbH (Munich); in South Pacific, contact Gammasonics Pty. Ltd. (Australia); in Japan, contact Toyo Medic Co. Ltd. (Tokyo).

² In U.S., contact Biodex Medical Systems (Shirley, New York); internationally, contact the U.S. office to find the dealer closest to you.

³ Flammability hazard. Do not use in or near operating centrifuges.

Storage

When the rotor is not in use, store it in a dry environment (not in the instrument) with the lid removed to allow air circulation so moisture will not collect in the tube cavities.

Returning a Rotor

Before returning a rotor or accessory for any reason, prior permission (a Returned Goods Authorization form) must be obtained from Beckman Coulter, Inc. This RGA form may be obtained from your local Beckman Coulter sales office. It should contain the following information:

- serial number,
- history of use (approximate frequency of use),
- reason for the return,
- original purchase order number, billing number, and shipping number, if possible,
- name and phone number of the person to be notified upon receipt of the rotor or accessory at the factory, and
- name and phone number of the person to be notified about repair costs, etc.

To protect our personnel, it is the customer's responsibility to ensure that the parts are free from pathogens and/or radioactivity. Sterilization and decontamination must be done before returning the parts. Smaller items (such as tubes, bottles, etc.) should be enclosed in a sealed plastic bag.

All parts must be accompanied by a note, plainly visible on the outside of the box or bag, stating that they are safe to handle and that they are not contaminated with pathogens or radioactivity. **Failure to attach** this notification will result in return or disposal of the items without review of the reported problem.

Use the address label printed on the RGA form when mailing the rotor and/or accessories.

Customers located outside the United States should contact their local Beckman Coulter office.

1-20 PN J-TB-035AT

Supply List

NOTE Publications referenced in this manual can be obtained at www.beckman.com, by calling Beckman Coulter at 1-800-742-2345 in the United States, or by contacting your local Beckman Coulter office.

See the *Beckman Coulter High Performance*, *High Speed*, *High Capacity Rotors*, *Tubes*, *and Accessories* catalog (PN BR-8102, available at www.beckman.com) for detailed information on ordering parts and supplies, or contact us. For your convenience, a partial list is given below.

Replacement Rotor Parts

JA-18 rotor assembly with magnets	
(rotors manufactured Jan. 2002 and later)	369679
JA-18 rotor assembly (rotors manufactured through Dec. 2001)	
Rotor lid assembly	364845
Rotor tie-down kit	364917

Other

Tubes, bottles, and accessories	see Table 1
Quick-Seal Cordless Tube Topper kit, 60 Hz	
Quick-Seal Cordless Tube Topper kit, 50 Hz (Europe)	
Quick-Seal Cordless Tube Topper kit, 50 Hz (Great Britain)	
Quick-Seal Cordless Tube Topper kit, 50 Hz (Australia)	
Quick-Seal Cordless Tube Topper kit, 50 Hz (Canada)	
Tube Topper rack, 12 place	
Spinkote lubricant (2 oz)	
Silicone vacuum grease (1 oz)	335148
Rotor Cleaning Kit	
Beckman Solution 555 (1 qt)	
Rotor cleaning brush	
Rotor removal tool and lifting handle	346965

PN J-TB-035AT 1-21

JA-18 Fixed Angle Rotor Supply List

1-22 PN J-TB-035AT

Beckman Coulter, Inc. J Series Rotor Warranty

Subject to the conditions specified below and the warranty clause of the Beckman Coulter, Inc., terms and conditions of sale in effect at the time of sale, Beckman Coulter, Inc. agrees to correct either by repair, or, at its election, by replacement, any defects of material or workmanship which develop within seven (7) years after delivery of a J series rotor to the original buyer by Beckman Coulter, Inc. or by an authorized representative, provided that investigation and factory inspection by Beckman Coulter, Inc. discloses that such defect developed under normal and proper use. Should a Beckman Coulter centrifuge be damaged due to a failure of a rotor covered by this warranty, Beckman Coulter will supply free of charge all centrifuge parts required for repair.

Replacement

Any product claimed to be defective must, if requested by Beckman Coulter, Inc., be returned to the factory, transportation charges prepaid, and will be returned to Buyer with the transportation charges collect unless the product is found to be defective, in which case Beckman Coulter, Inc. will pay all transportation charges.

A defective rotor will be replaced by Beckman Coulter, Inc. at its then current list price less a credit based upon the age of the rotor (years since date of purchase). The Buyer shall not receive credit until the claimed defective rotor is returned to Beckman Coulter's Indianapolis, Indiana facility or delivered to a Beckman Field Service representative.

The replacement price (cost to Buyer) for the respective rotor shall be calculated as follows:

Replacement price = Current rotor list price $\times \frac{\text{years}}{7}$

Conditions

- 1. Except as otherwise specifically provided herein, this warranty covers the rotor only and Beckman Coulter, Inc. shall not be liable for damage to accessories or ancillary supplies including but not limited to (i) tubes, (ii) tube caps, (iii) tube adapters, or (iv) tube contents.
- **2.** This warranty is void if the rotor has been subjected to customer misuse such as operation or maintenance contrary to the instructions in the Beckman Coulter rotor or centrifuge manual.
- **3.** This warranty is void if the rotor is operated with a rotor drive unit or in a centrifuge unmatched to the rotor characteristics, or is operated in a Beckman Coulter centrifuge that has been improperly disassembled, repaired, or modified.

Disclaimer

IT IS EXPRESSLY AGREED THAT THE ABOVE WARRANTY SHALL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND OF THE WARRANTY OF MERCHANTABILITY AND THAT BECKMAN COULTER, INC. SHALL HAVE NO LIABILITY FOR SPECIAL OR CONSEQUENTIAL DAMAGES OF ANY KIND WHATSOEVER ARISING OUT OF THE MANUFACTURE, USE, SALE, HANDLING, REPAIR, MAINTENANCE, OR REPLACEMENT OF THE PRODUCT.

PN J-TB-035AT Warranty-1

Beckman Coulter, Inc. J Series Rotor Warranty

Warranty-2 PN J-TB-035AT

Related Documents

Rotors and Tubes for Beckman Coulter J2, J6, and Avanti J Series Centrifuges

PN JR-IM-10

- Rotors
- Tubes, Bottles, and Accessories
- Using Tubes and Accessories
- Using Fixed-Angle Rotors
- Using Swinging-Bucket Rotors
- Using Vertical-Tube and Rack Type Rotors
- Care and Maintenance
- Chemical Resistances
- Temperature Compensation Tables
- Gradient Materials
- Blood Component Separation

Use and Care of Centrifuge Tubes and Bottles

PN IN-192

Chemical Resistances for Beckman Coulter Centrifugation Products

PN IN-175

High Performance, High Speed, High Capacity Rotors, Tubes, & Accessories

PN BR-1802

www.beckman.com

